

Fast Analysis of Non-Traditional Gasoline Additives with Gas Chromatography-Vacuum Ultraviolet Spectroscopy


Ryan Schonert, Dan Wispinski, Jack Cochran






Gasoline Additives



- Organic compounds which adjust fuel properties
 - Refining process, store-bought
- Major fuel blending components: octane boosters, pollution preventers
 - Blended at volume % level

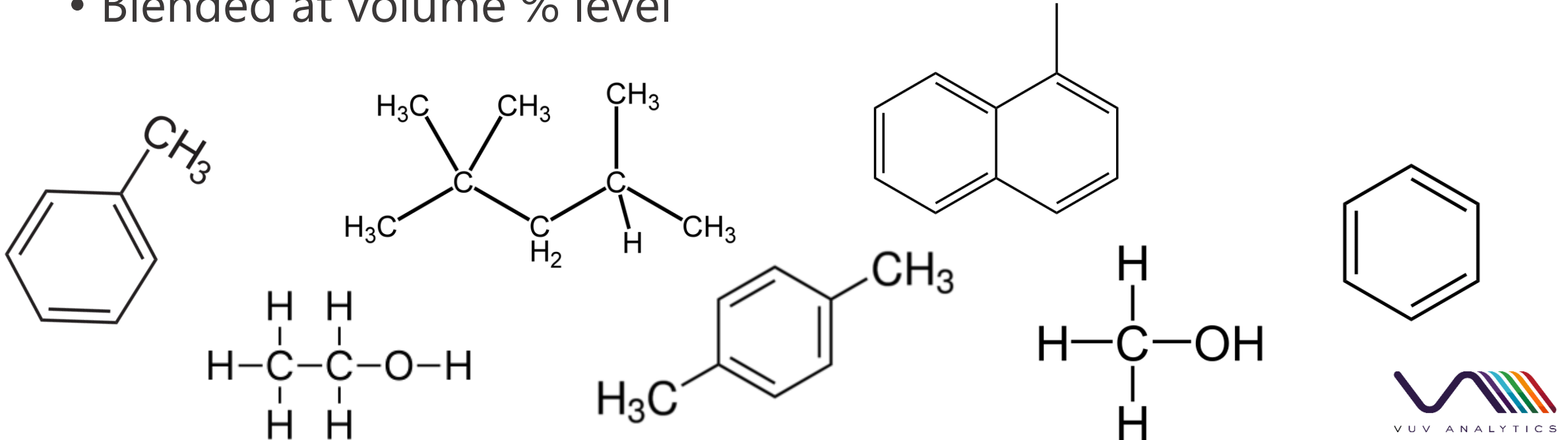

Gasoline Additives



- Organic compounds which adjust fuel properties
 - Refining process, store-bought
- Major fuel blending components: octane boosters, pollution preventers
 - Blended at volume % level


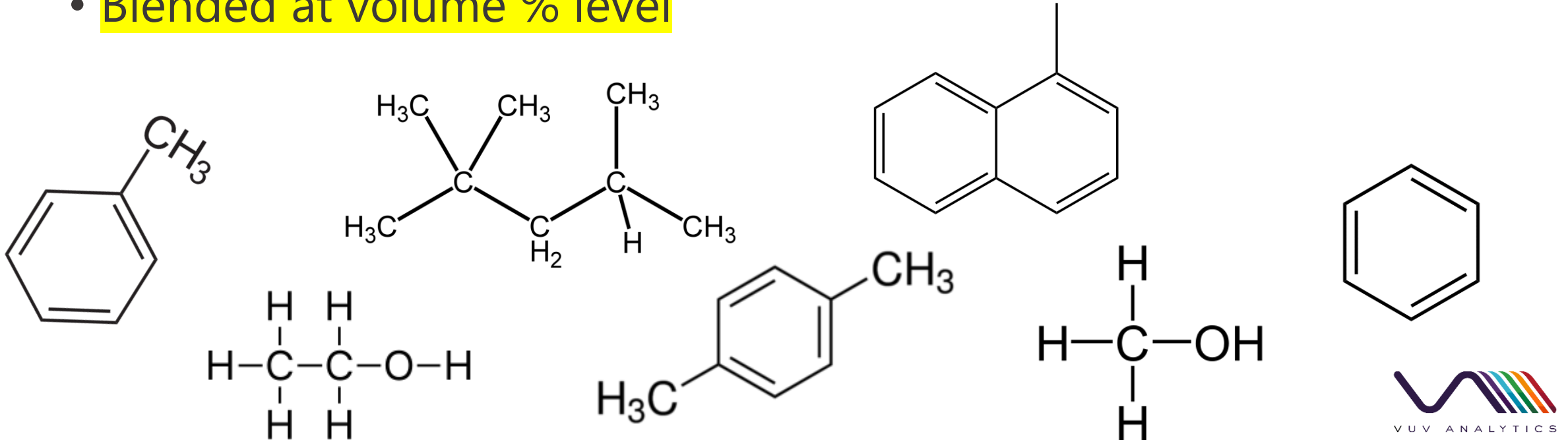


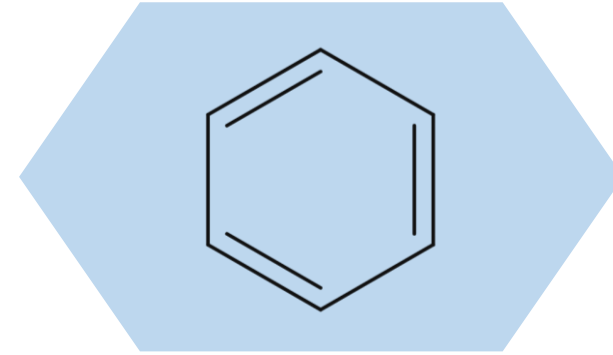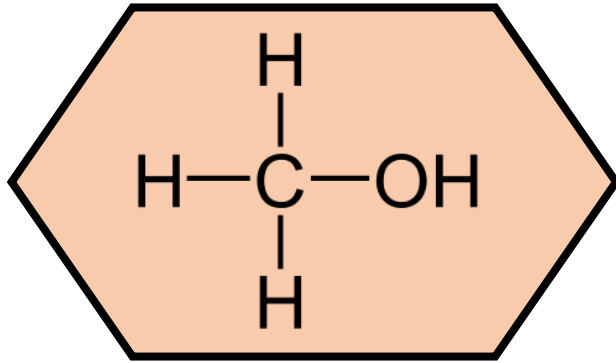

Methanol

- Cheap and efficient octane booster
- Emissions: high formaldehyde concentrations

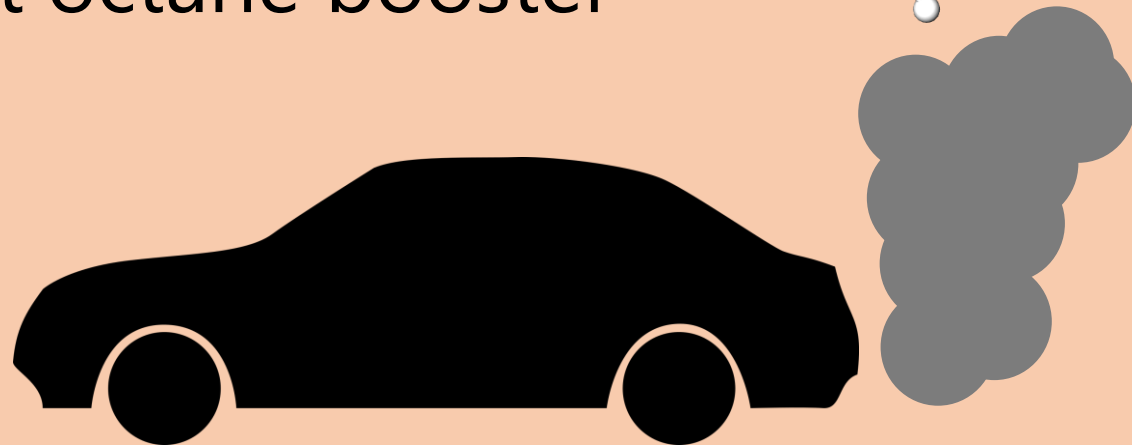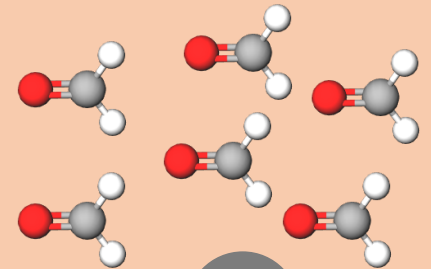

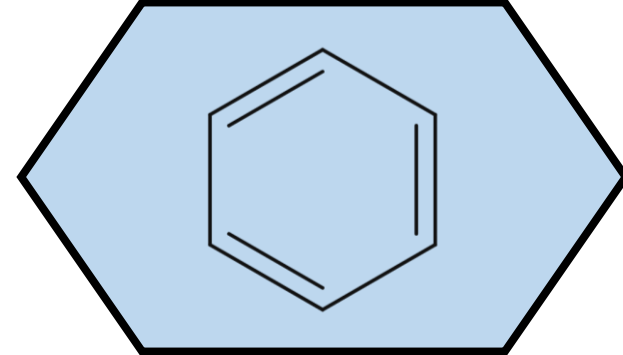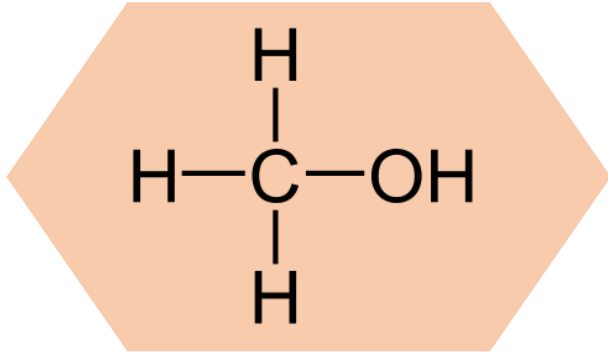

Benzene

- Important BTEX octane-boosting component
- Group 1 carcinogen



Non-Traditional Gasoline Additives (NTGAs)


Desired Effect



Environmentally Safe




Renewable

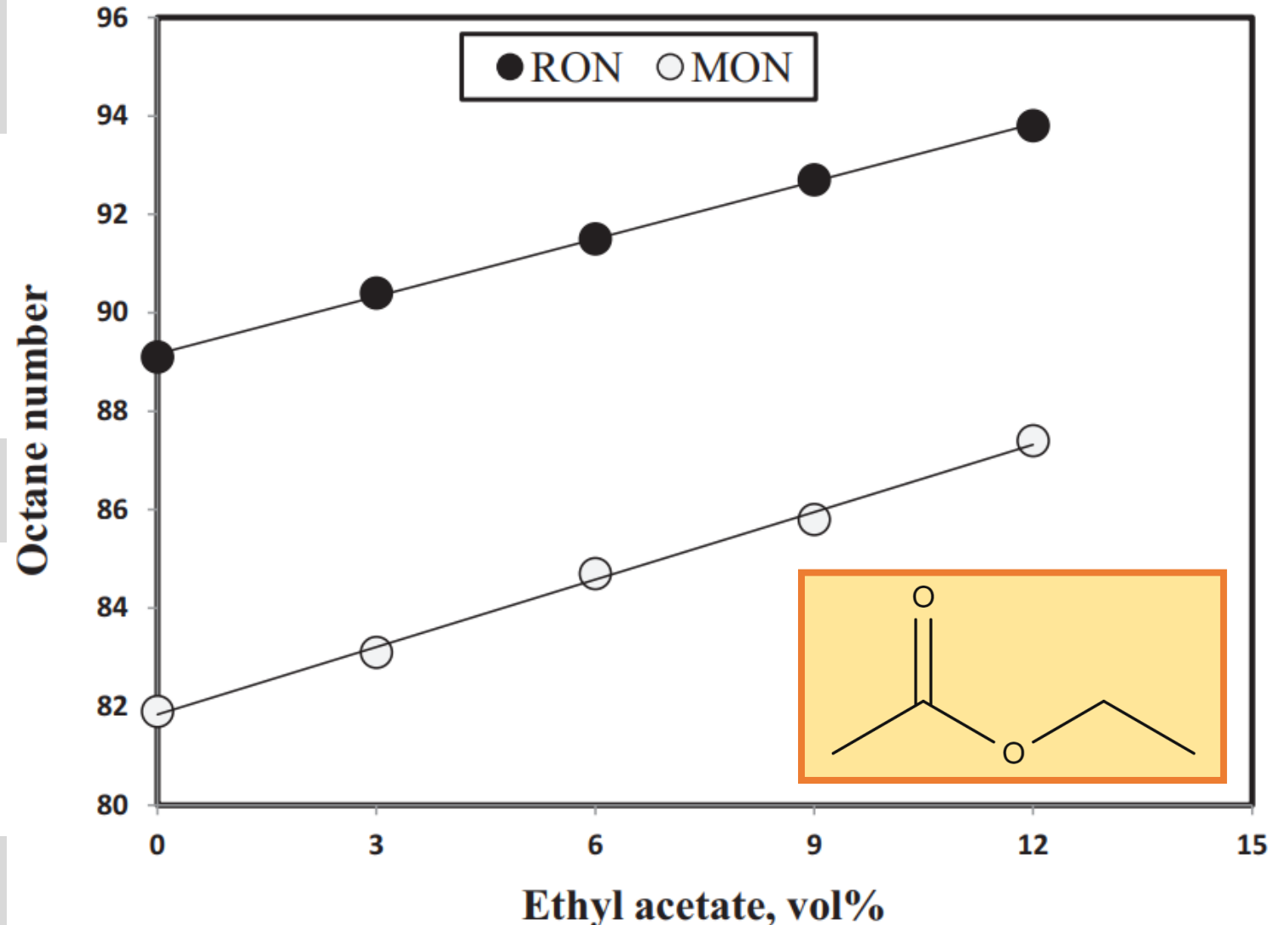



Home > Vol 5, No 2 > Irías-Mata

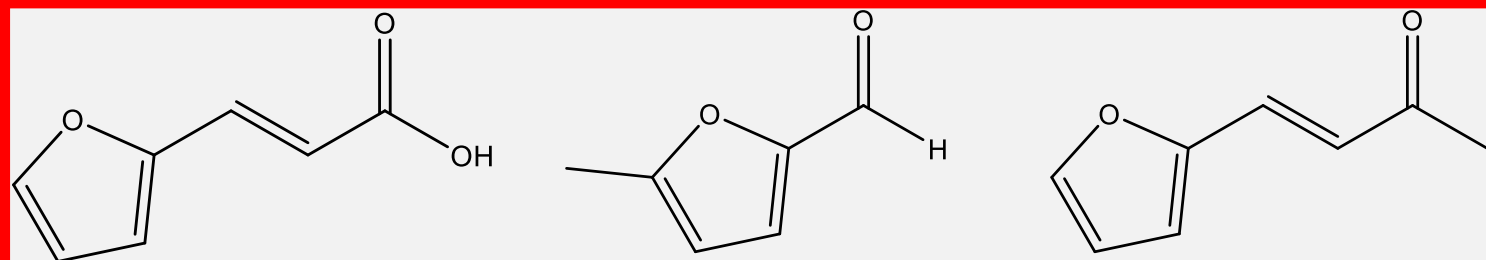

Pineapple-stover derived furan compounds as gasoline oxygenate additive

Andrea P. Irías-Mata, Giselle Lutz

Abstract

Furan compounds have properties such as oxygenate additive to enhance octane number of gasoline. The procedure was a furan synthesis through an acidic hydrolysis of the polysaccharide materials from pineapple plantation residues. The products obtained were a complex mixture of mostly 3-(2-furanyl)-2-propenoic acid, 4-(2-furanyl)-3-butene-2-one and 5-methyl-furfural. Thermodynamic and rheological properties of the mixture *in toto* were measured, as well as its oxygenating capability. The results showed a two units enhancer oxygenate additive for naphta, potentially safe for transport and handling, presenting the following characteristics: specific gravity $1,22559 \pm 0,00002$, kinematic viscosity $0,0127 \pm 0,0001$ Stokes, enthalpy of vaporization $39,1 \pm 0,1 \text{ kJ mol}^{-1}$, isothermal compressibility $(1,0 \pm 0,2) 10^{-9} \text{ Pa}^{-1}$, rate of evaporation $0,03 \pm 0,02 \text{ g s}^{-1} \text{ m}^{-2}$ and Hildebrand solubility parameter $18,0 \pm 0,1 (\text{J cm}^{-3})^{1/2}$.

KEY WORDS

Pineapple crown, pineapple stubble, research octane number, oxygenate additive, furan compound, methyl terbutyl ether.

Selection Criteria and Screening of Potential Biomass-Derived Streams as Fuel Blendstocks for Advanced Spark-Ignition Engines

Robert L. McCormick, Gina Fioroni, Lisa Fouts, and Earl Christensen
National Renewable Energy Laboratory

Janet Yanowitz
Ecoengineering Inc.

Evgueni Polikarpov, Karl Albrecht, and Daniel J. Gaspar
Pacific Northwest National Laboratory

John Gladden and Anthe George
Sandia National Laboratory

***Biomass-derived
compounds that
have potential as
fuel blendstocks
(partial list)***

Name	Formula	CAS#	RON (D2699)	MON (D2700)	HOV, kJ/kg ^a	LHV, MJ/kg	LHV, MJ/L (at 20°C)	Density, g/L (at 20°C)	Water Solubility, mg/L
<i>Alcohols</i>									
Methanol	CH ₃ O	67-56-1	109	89	1173.5	20.09	15.8	786.4	miscible
Ethanol	C ₂ H ₆ O	64-17-5	109	90	918.2	25.54	20.2	789.3	miscible
n-Propanol ^c	C ₃ H ₈ O	71-23-8	104	89	788.7	30.8	24.7	811.0	miscible
2-propanol ^c	C ₃ H ₈ O	67-63-0	109	97	743.8	30.7	24.1	781.2	miscible
1-Butanol	C ₄ H ₁₀ O	71-36-3	98	85	708.3	33.1	26.8	805.7	68,000
2-butanol	C ₄ H ₁₀ O	78-92-2	107	93	670.5	33.1	26.7	802.6	181,000
2-methylpropan-1-ol (isobutanol)	C ₄ H ₁₀ O	78-83-1	105	90	685.4	33.2	26.6	797.6	85,000
2-methyl-1-butanol	C ₅ H ₁₂ O	137-32-6	101	88.3	613.7	34.69	28.3	814.8	29,700
2-pentanol	C ₅ H ₁₂ O	6032-29-7	99.4	90.8	608.1	34.58	27.8	805.2	44,600
<i>Alkanes</i>									
Triptane (2,2,3-trimethylbutane) ^c	C ₇ H ₁₆	464-06-2	112	101	319.3	44.4	30.8	694.5	29
<i>Alkenes</i>									
Diisobutylene (a mixture of three parts 2,4,4-trimethyl-1-pentene and one part 2,4,4-trimethyl-2-pentene)	C ₈ H ₁₆	25167-70-8	106	86.5	318.2	44.27	31.66	715.1	4
<i>Esters</i>									
Methylacetate	C ₃ H ₆ O ₂	79-20-9	>120	>120	438.7	19.3	17.9	927.1	243,000
Methylbutanoate ^c	C ₅ H ₁₀ O ₂	623-42-7	107.2	105	401.4	26.07	23.25	892.2	15,000
Methylisobutanoate (2-methylpropanoate) ^c	C ₅ H ₁₀ O ₂	547-63-7	103.6	104.7	365.2	28.38	25.04	882.4	9,260
Methylpentanoate ^c	C ₆ H ₁₂ O ₂	624-24-8	103.4	101.1	379.6	28.25	24.99	884.8	2,200
Methyl-2-methylbutanoate ^c	C ₆ H ₁₂ O ₂	868-57-5	110.5	99.1	338.3 ^b	28.08	24.59	875.7	3,200
Ethylacetate	C ₄ H ₈ O ₂	141-78-6	118	>120	399.5	23.79	21.34	894.6	80,000
Ethylbutanoate	C ₆ H ₁₂ O ₂	105-54-4	115.4	106	369.3	28.64	25.17	874.3	4,900
Ethylisobutyrate (2-methylpropanoate) ^c	C ₆ H ₁₂ O ₂	97-62-1	110.3	109.5	342.6	30.44	26.46	869.3	3,200
1-Methylethylacetate (isopropylacetate) ^c	C ₅ H ₁₀ O ₂	108-21-4	>120	--	363.3	25.77	22.47	871.8	29,000
Butylacetate	C ₆ H ₁₂ O ₂	123-86-4	100.8	100	369.3	27.43	24.03	876.0	8,400
Isobutylacetate (2-methylpropylacetate) ^c	C ₆ H ₁₂ O ₂	110-19-0	108.7	112.3	319.4	28.01	24.4	871.2	6,300
Isoamylacetate (3-methylbutylacetate) ^c	C ₇ H ₁₄ O ₂	123-92-2	100.6	97.2	356.4	32.15	27.84	866.1	2,000
<i>Ethers</i>									
Methoxybenzene (anisole)	C ₇ H ₈ O	100-66-3	103	92	428.2	33.5	33.3	989.2	1,040
<i>Furans</i>									
2-Methylfuran/2,5-dimethylfuran mixture (40/60 by weight)	--	--	102	87		32.2	29.1	903	2,240
<i>Ketones</i>									
2-Butanone (methyl ethyl ketone)	C ₄ H ₈ O	78-93-3	111	105.5	481.2	31.36	25.08	799.8	223,000
2-Pentanone	C ₅ H ₁₀ O	107-87-9	105.7	103	445.8	33.37	26.75	801.5	43,000
3-Pentanone ^c	C ₅ H ₁₀ O	96-22-0	106.8	--	448.2	33.43	27.08	810.1	45,900
3-Methyl-2-butanone ^c		563-80-4	108.9	102.2	428.4	33.34	26.68	815.0	5,237
Cyclopentanone	C ₅ H ₈ O	120-92-3	101	89.4	504.0	31.99	30.2	944.1	60,000
3-Hexanone ^c	C ₆ H ₁₂ O	589-38-8	101.9	93.5	417.3	34.83	28.61	821.6	14,700
4-Methyl-2-pentanone (methylisobutyl ketone) ^c	C ₆ H ₁₂ O	108-10-1	105.7	105.5	417.3	34.21	27.27	797.1	19,000
2,4-dimethyl-3-pentanone	C ₇ H ₁₄ O	565-80-0	99	92.5	363.4	35.84	28.81	803.9	5,700
Ketone mixture (42.5 wt% 2-pentanone, 11.4 wt% methyl-isobutyl ketone, 30.3 wt% 4-heptanone, 15.8 wt% 2-heptanone)	--	--	99.4	99.6	424 ^b	34.61	27.99	808.7	--

Selection Criteria and Screening of Potential Biomass-Derived Streams as Fuel Blendstocks for Advanced Spark-Ignition Engines

Robert L. McCormick, Gina Fioroni, Lisa Fouts, and Earl Christensen

Alcohols

Alkanes

Alkenes

Esters

Ethers

Furans

Ketones

Name	Formula	CAS#	RON (D2699)	MON (D2700)	HOV, kJ/kg ^a	LHV, MJ/kg	LHV, MJ/L (at 20°C)	Density, g/L (at 20°C)	Water Solubility, mg/L
<i>Alcohols</i>									
Methanol	CH ₄ O	67-56-1	109	89	1173.5	20.09	15.8	786.4	miscible
Ethanol	C ₂ H ₆ O	64-17-5	109	90	918.2	25.54	20.2	789.3	miscible
n-Propanol ^c	C ₃ H ₈ O	71-23-8	104	89	788.7	30.8	24.7	811.0	miscible
2-propanol ^c	C ₃ H ₈ O	67-63-0	109	97	743.8	30.7	24.1	781.2	miscible
1-Butanol	C ₄ H ₁₀ O	71-36-3	98	85	708.3	33.1	26.8	805.7	68,000
2-butanol	C ₄ H ₁₀ O	78-92-2	107	93	670.5	33.1	26.7	802.6	181,000
2-methylpropan-1-ol (isobutanol)	C ₄ H ₁₀ O	78-83-1	105	90	685.4	33.2	26.6	797.6	85,000
2-methyl-1-butanol	C ₅ H ₁₂ O	137-32-6	101	88.3	613.7	34.69	28.3	814.8	29,700
2-pentanol	C ₅ H ₁₂ O	6032-29-7	99.4	90.8	608.1	34.58	27.8	805.2	44,600
<i>Alkanes</i>									
Triptane (2,2,3-trimethylbutane) ^c	C ₇ H ₁₆	464-06-2	112	101	319.3	44.4	30.8	694.5	29
<i>Alkenes</i>									
Diisobutylene (a mixture of three parts 2,4,4-trimethyl-1-pentene and one part 2,4,4-trimethyl-2-pentene)	C ₈ H ₁₆	25167-70-8	106	86.5	318.2	44.27	31.66	715.1	4
<i>Esters</i>									
Methylacetate	C ₃ H ₆ O ₂	79-20-9	>120	>120	438.7	19.3	17.9	927.1	243,000
Methylbutanoate ^c	C ₅ H ₁₀ O ₂	623-42-7	107.2	105	401.4	26.07	23.25	892.2	15,000
Methylisobutanoate (2-methylpropanoate) ^c	C ₅ H ₁₀ O ₂	547-63-7	103.6	104.7	365.2	28.38	25.04	882.4	9,260
Methylpentanoate ^c	C ₆ H ₁₂ O ₂	624-24-8	103.4	101.1	379.6	28.25	24.99	884.8	2,200
Methyl-2-methylbutanoate ^c	C ₆ H ₁₂ O ₂	868-57-5	110.5	99.1	338.3 ^b	28.08	24.59	875.7	3,200
Ethylacetate	C ₄ H ₈ O ₂	141-78-6	118	>120	399.5	23.79	21.34	894.6	80,000
Ethylbutanoate	C ₆ H ₁₂ O ₂	105-54-4	115.4	106	369.3	28.64	25.17	874.3	4,900
Ethylisobutyrate (2-methylpropanoate) ^c	C ₆ H ₁₂ O ₂	97-62-1	110.3	109.5	342.6	30.44	26.46	869.3	3,200
1-Methylethylacetate (isopropylacetate) ^c	C ₅ H ₁₀ O ₂	108-21-4	>120	--	363.3	25.77	22.47	871.8	29,000
Butylacetate	C ₆ H ₁₂ O ₂	123-86-4	100.8	100	369.3	27.43	24.03	876.0	8,400
Isobutylacetate (2-methylpropylacetate) ^c	C ₆ H ₁₂ O ₂	110-19-0	108.7	112.3	319.4	28.01	24.4	871.2	6,300
Isoamylacetate (3-methylbutylacetate) ^c	C ₇ H ₁₄ O ₂	123-92-2	100.6	97.2	356.4	32.15	27.84	866.1	2,000
<i>Ethers</i>									
Methoxybenzene (anisole)	C ₇ H ₈ O	100-66-3	103	92	428.2	33.5	33.3	989.2	1,040
<i>Furans</i>									
2-Methylfuran/2,5-dimethylfuran mixture (40/60 by weight)	--	--	102	87		32.2	29.1	903	2,240
<i>Ketones</i>									
2-Butanone (methyl ethyl ketone)	C ₄ H ₈ O	78-93-3	111	105.5	481.2	31.36	25.08	799.8	223,000
2-Pentanone	C ₅ H ₁₀ O	107-87-9	105.7	103	445.8	33.37	26.75	801.5	43,000
3-Pentanone ^c	C ₅ H ₁₀ O	96-22-0	106.8	--	448.2	33.43	27.08	810.1	45,900
3-Methyl-2-butanone ^c		563-80-4	108.9	102.2	428.4	33.34	26.68	815.0	5,237
Cyclopentanone	C ₅ H ₈ O	120-92-3	101	89.4	504.0	31.99	30.2	944.1	60,000
3-Hexanone ^c	C ₆ H ₁₂ O	589-38-8	101.9	93.5	417.3	34.83	28.61	821.6	14,700
4-Methyl-2-pentanone (methylisobutyl ketone) ^c	C ₆ H ₁₂ O	108-10-1	105.7	105.5	417.3	34.21	27.27	797.1	19,000
2,4-dimethyl-3-pentanone	C ₇ H ₁₄ O	565-80-0	99	92.5	363.4	35.84	28.81	803.9	5,700
Ketone mixture (42.5 wt% 2-pentanone, 11.4 wt% methyl-isobutyl ketone, 30.3 wt% 4-heptanone, 15.8 wt% 2-heptanone)	--	--	99.4	99.6	424 ^b	34.61	27.99	808.7	--

Harmful NTGAs

Methyl Acetate

Acetone

N-Methylaniline

Methylal

Secondary-Butyl Acetate

“Considered to be a cheap source of octane, these chemicals could cause engine problems...[t]hey are also harmful to human health and to the environment...”

Harmful NTGAs


Methyl Acetate

N-Methylaniline

Secondary-Butyl Acetate

Acetone

Methylal

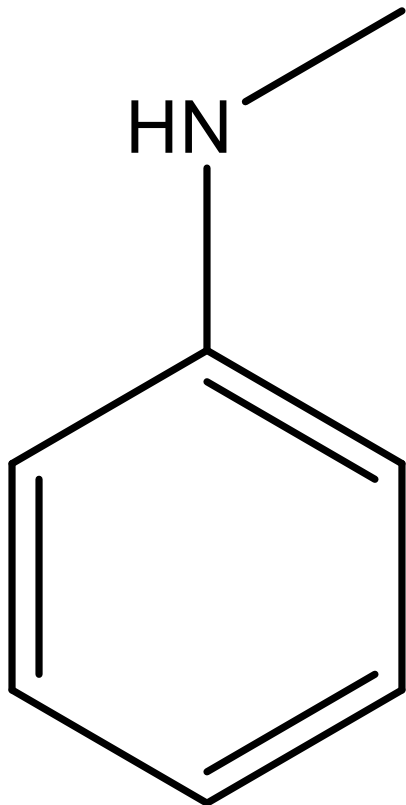

- **Octane booster**
- **Causes gum formation** – carbon deposit in engine parts
- **Swells rubber seals** – may cause oil leaks


Gasoline Analysis – DHA, ASTM D6730

- *2+ hour run time*
- *Identification by retention time only*
- ***Not equipped to analyze novel compounds***



Designation: D6730 – 01 (Reapproved 2016)

**Standard Test Method for
Determination of Individual Components in Spark Ignition
Engine Fuels by 100–Metre Capillary (with Precolumn) High-
Resolution Gas Chromatography¹**

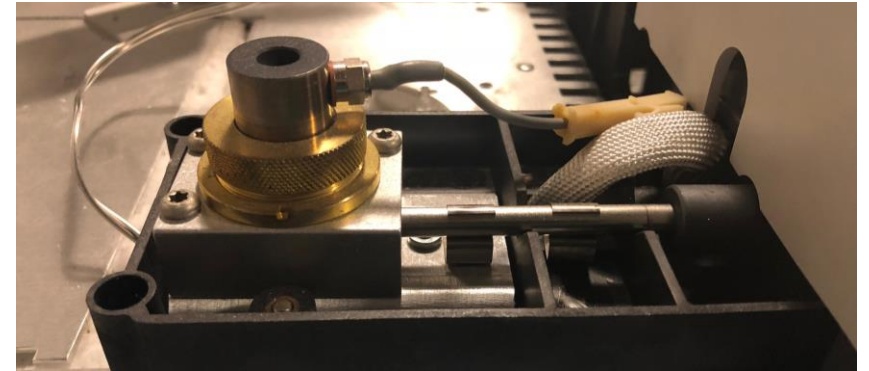


Gasoline Analysis – DHA, ASTM D6730

- 2+ hour run time
- Identification by retention time only
- **Not equipped to analyze novel compounds**




Designation: D6730 – 01 (Reapproved 2006)


Standard Test Method for
Determination of Individual Components in Spark Ignition
Engine Fuels by 100–Metre Capillary (with Precolumn) High-
Resolution Gas Chromatography¹

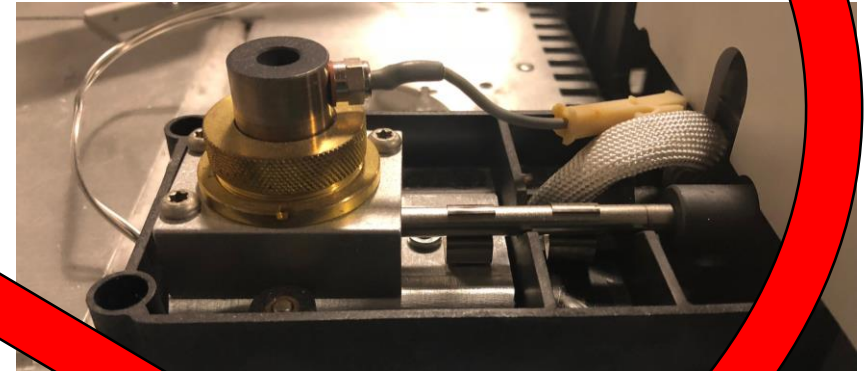


How Does VUV Spectroscopy Work?

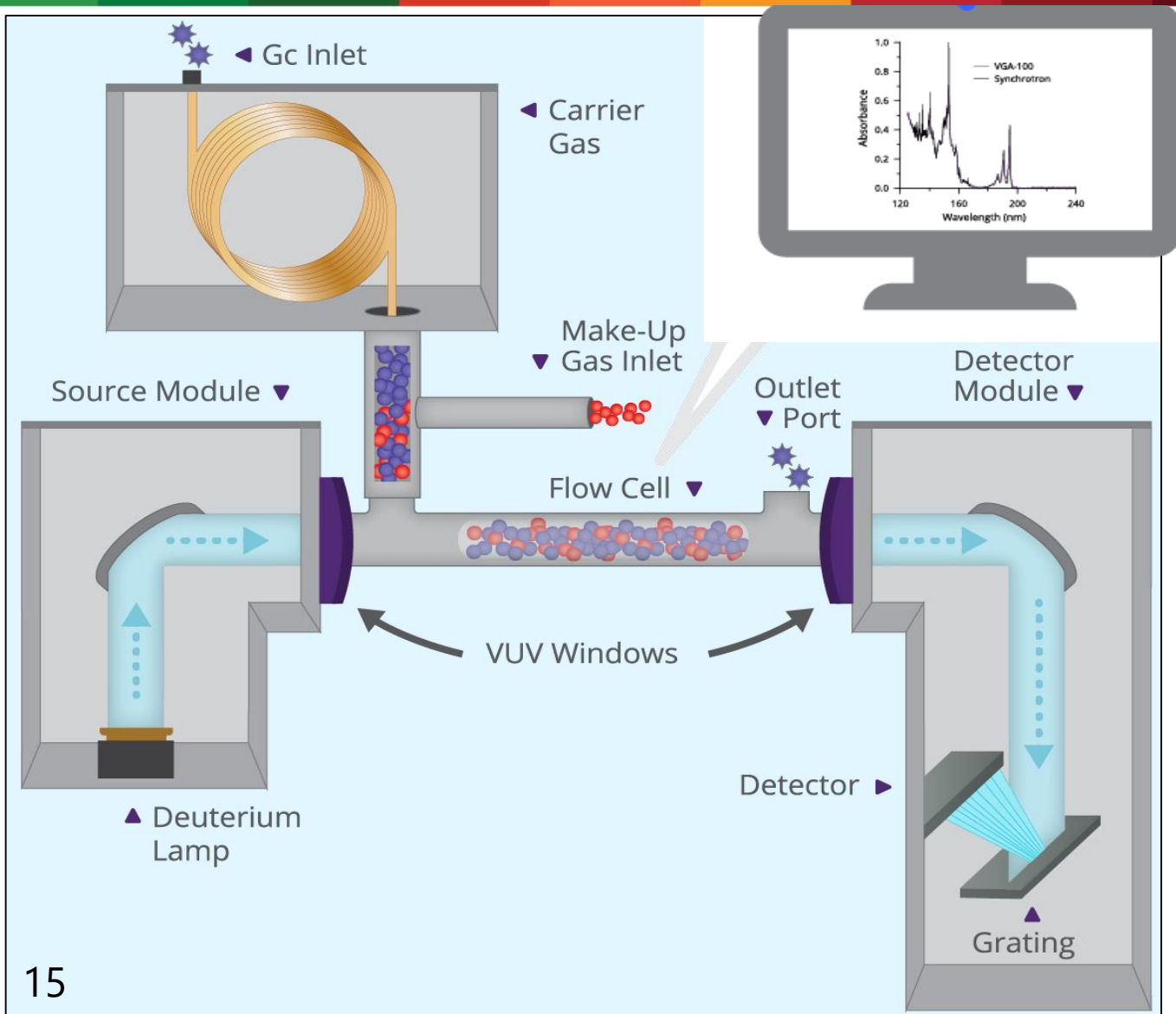


VUV Analytics VGA-100
125 to 240nm
1 to 75 spectra/sec




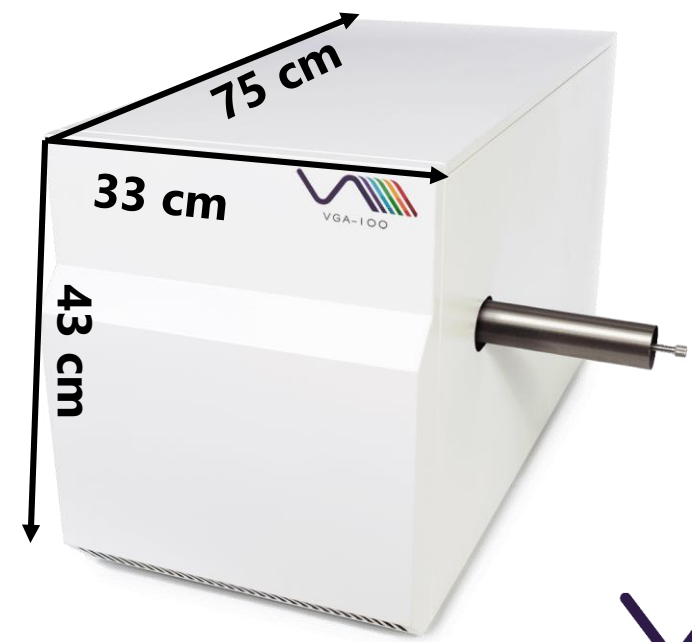



How Does VUV Spectroscopy Work?

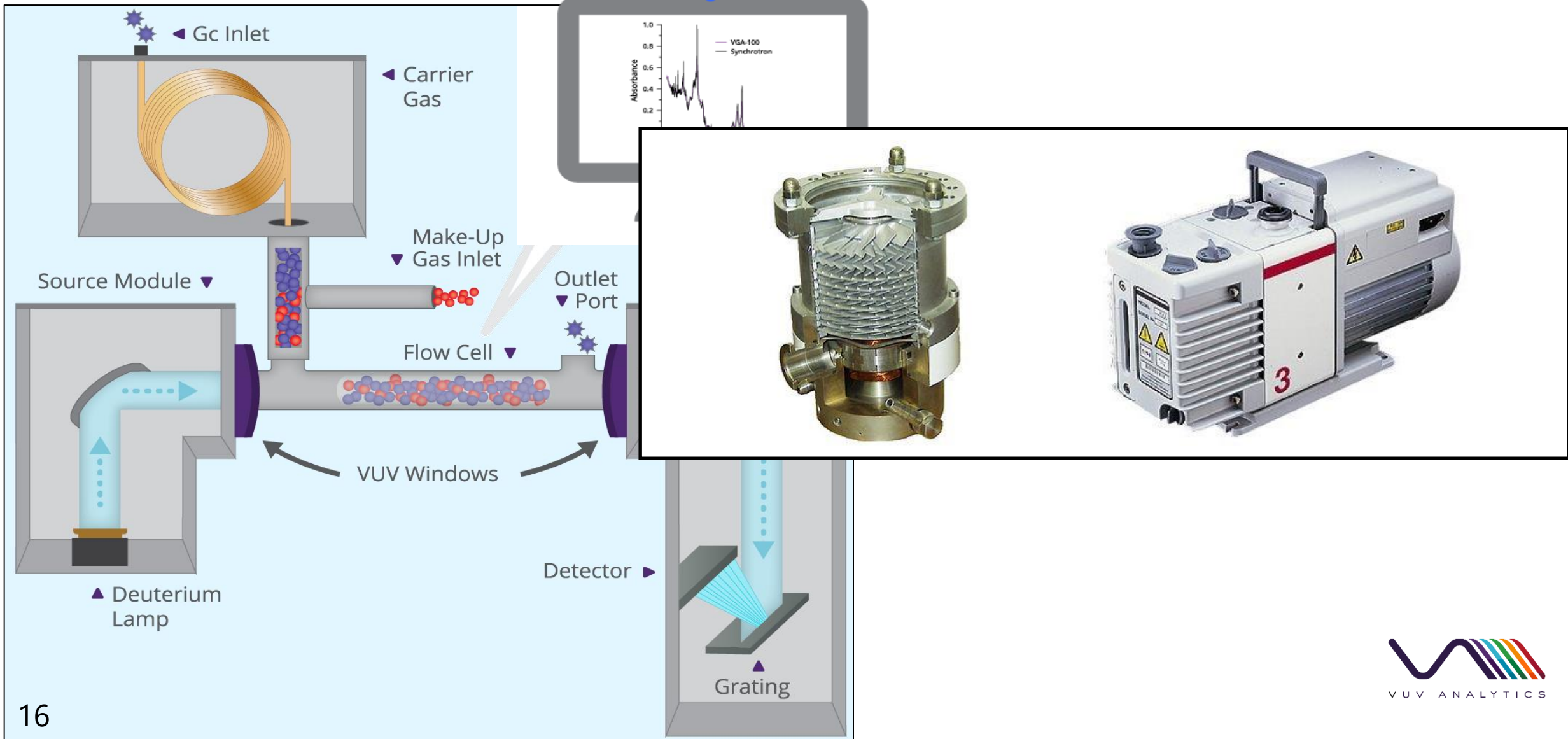


How Does VUV Spectroscopy Work?

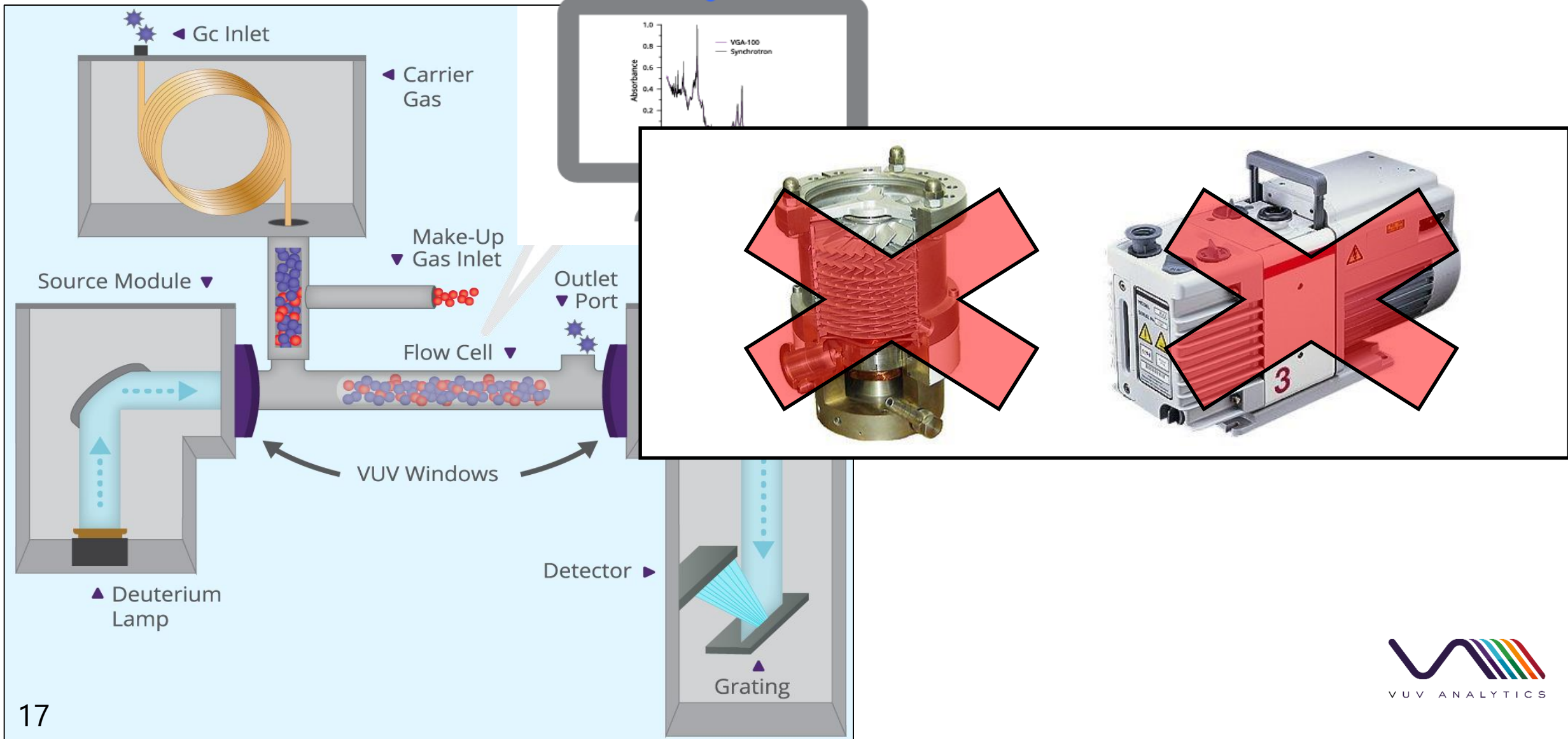

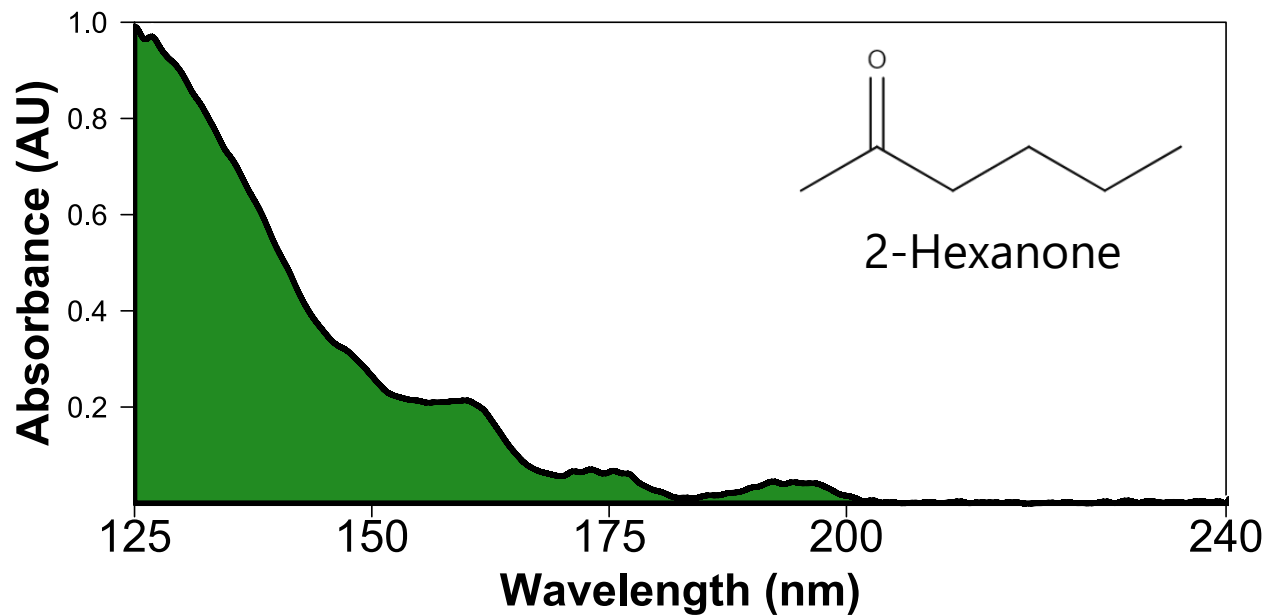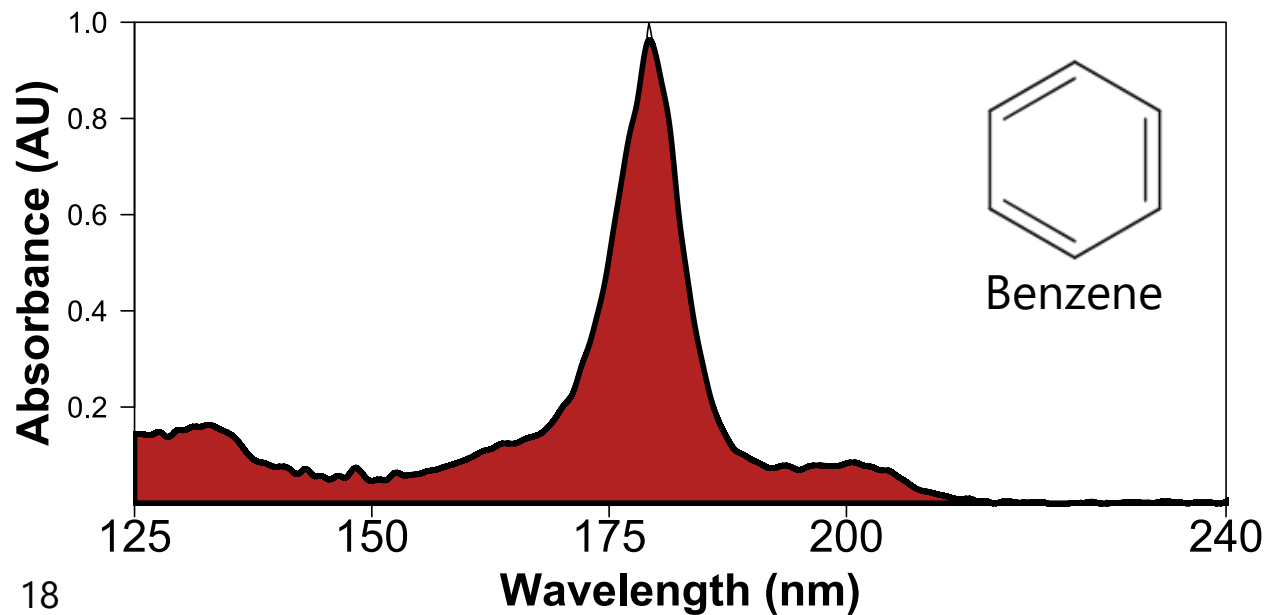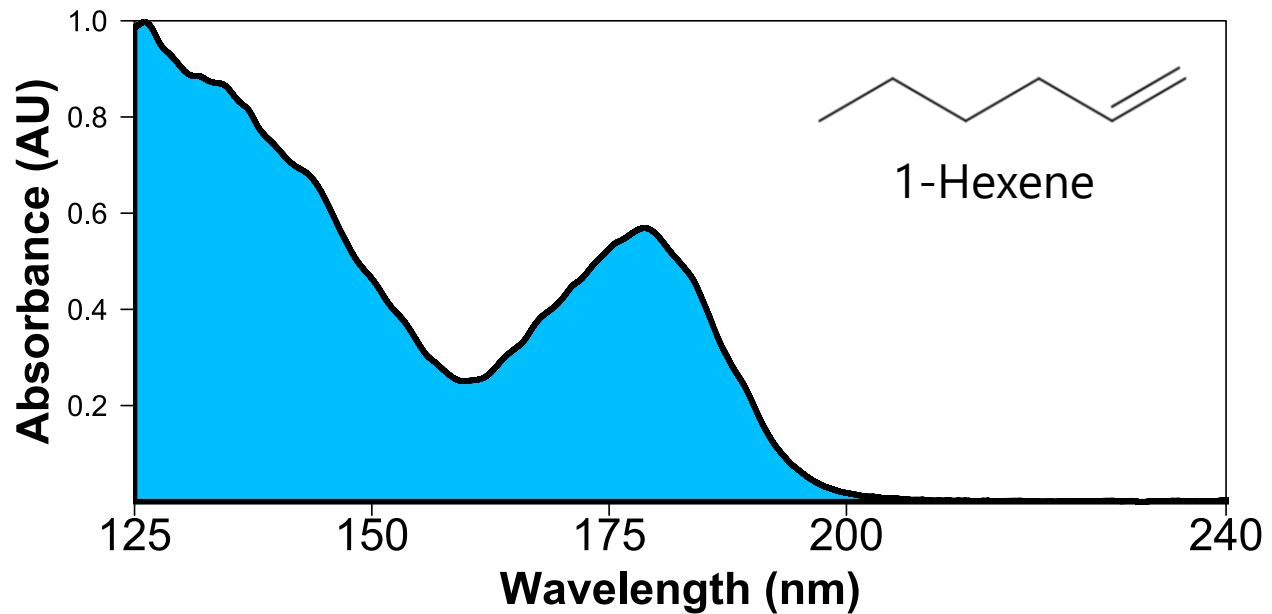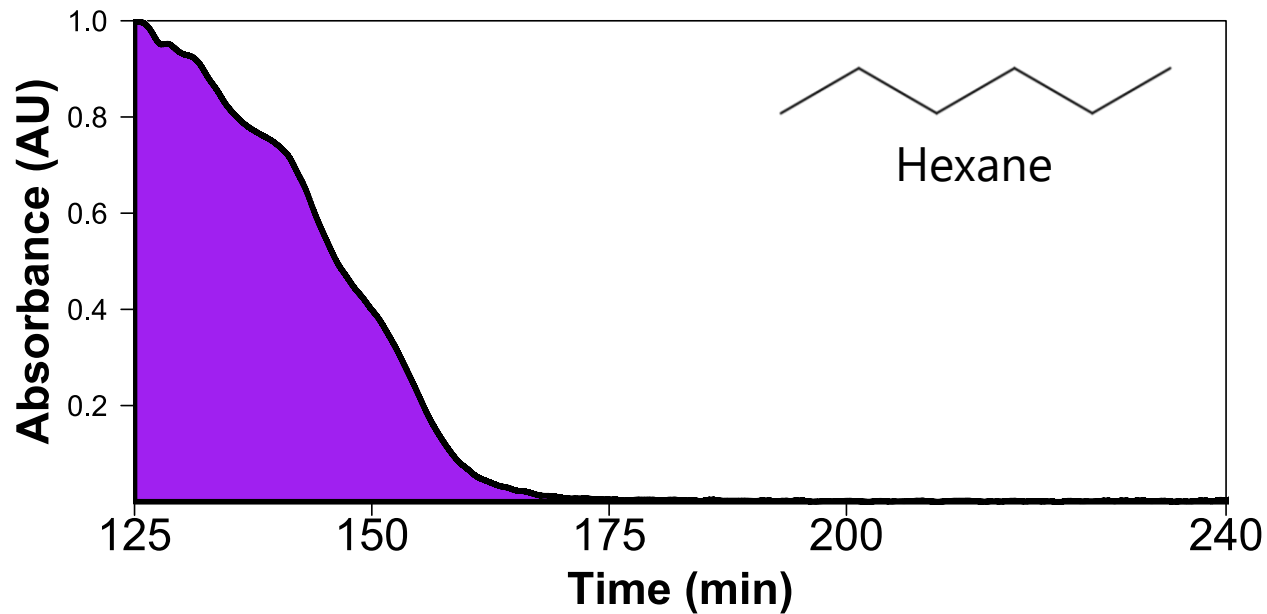


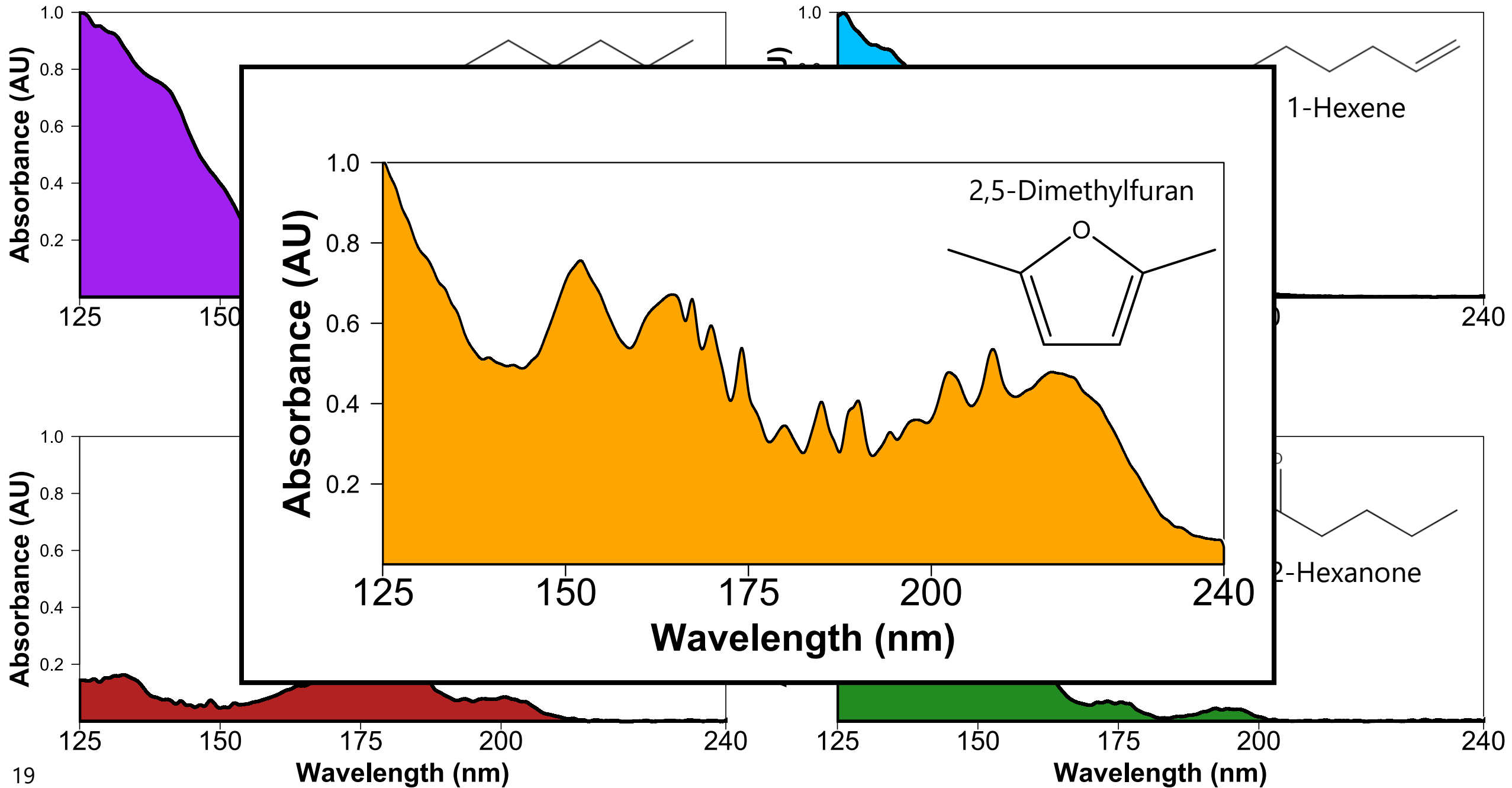


Coelution of Iso-pentane and Acetone

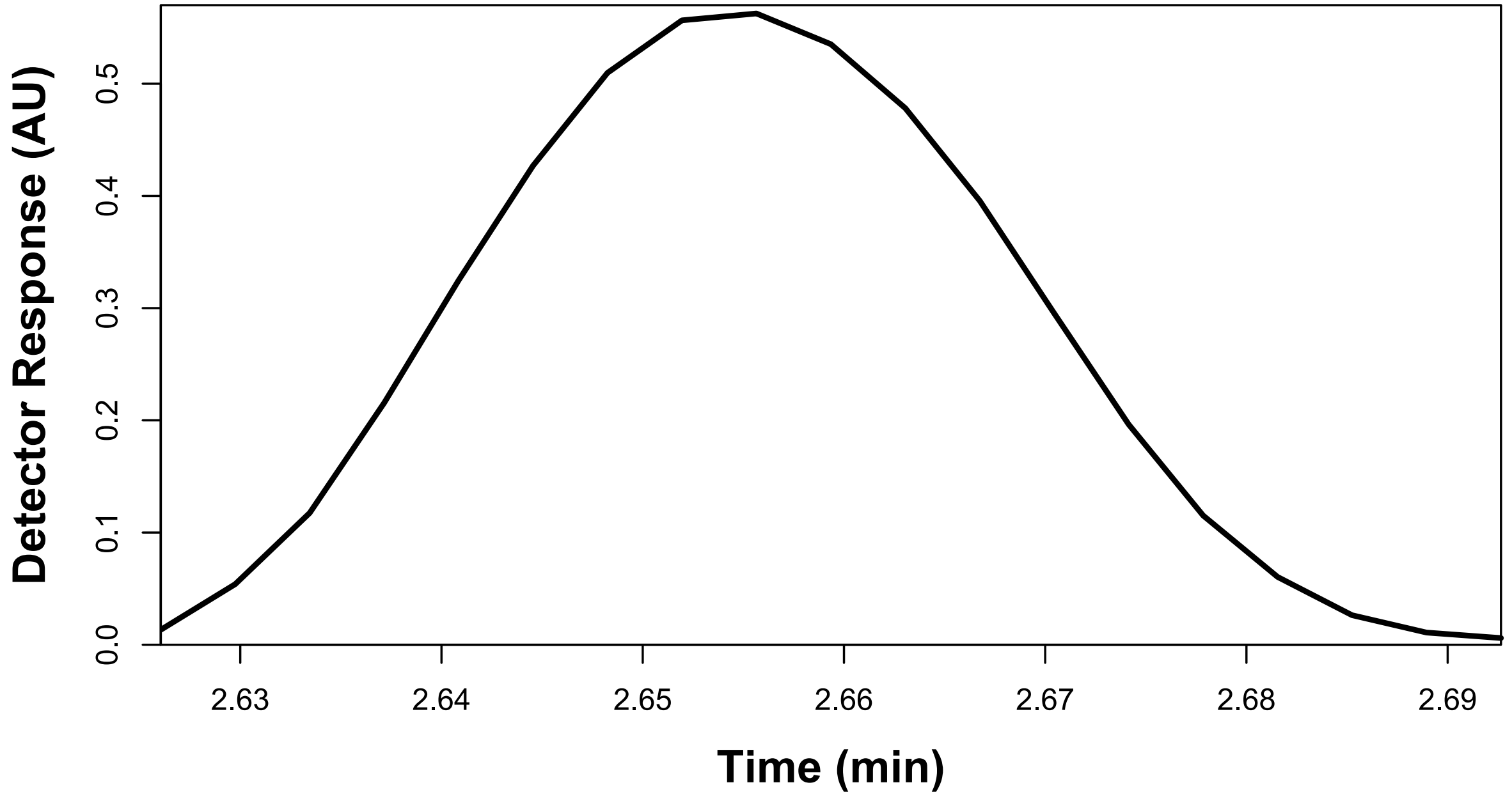


Coelution of Iso-pentane and Acetone

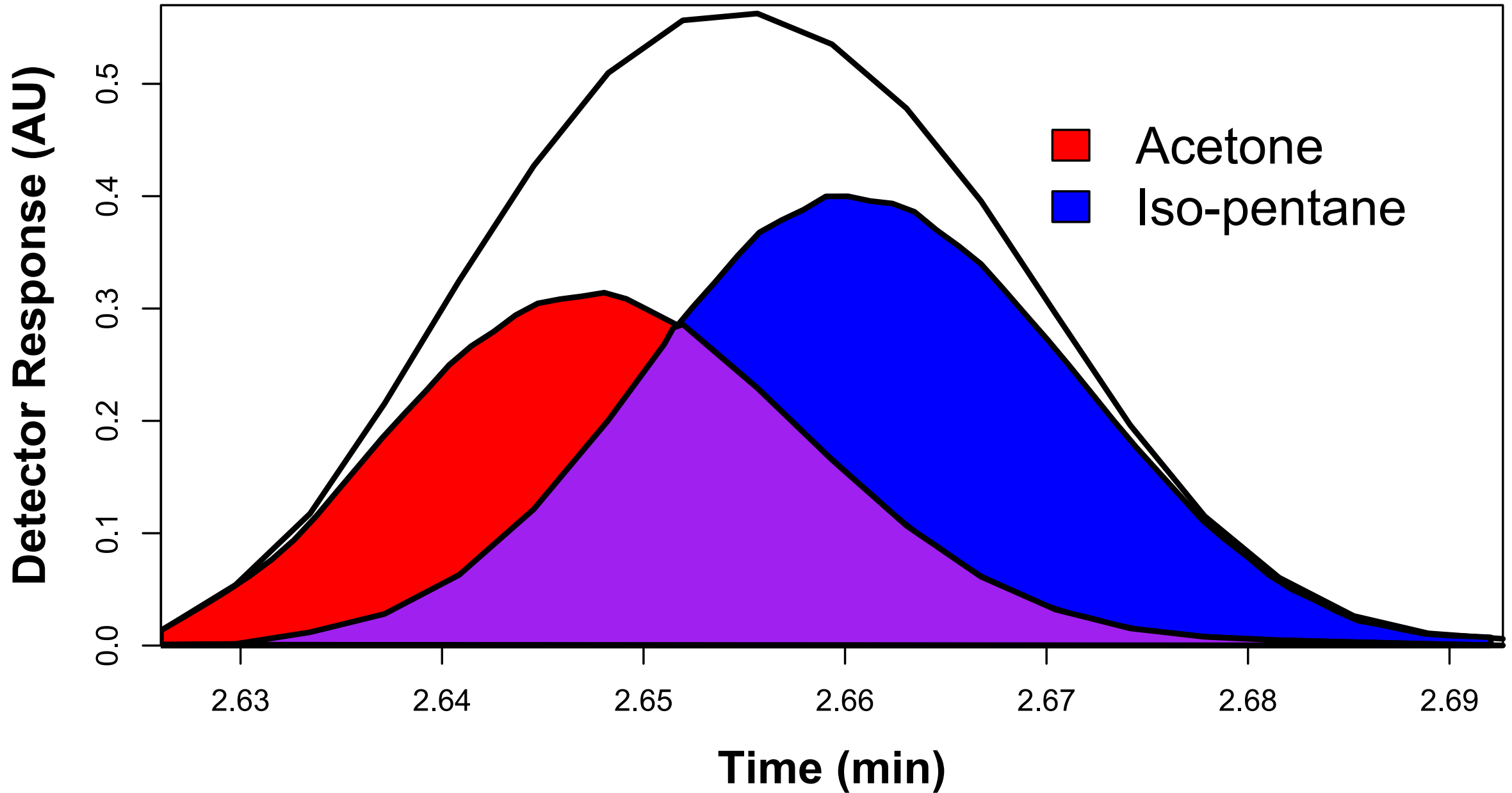


Coelution of Iso-pentane and Acetone

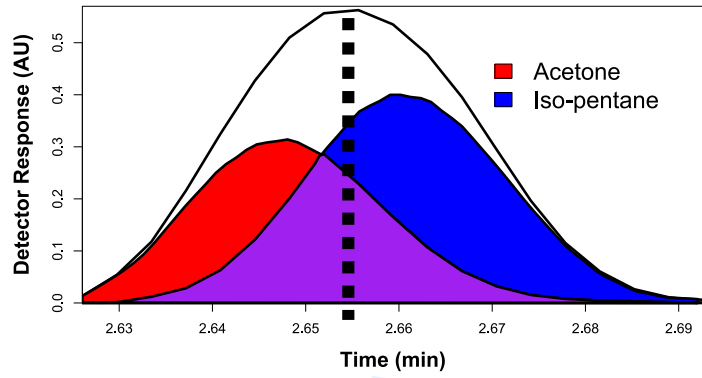


Spectral Absorbance of Acetone and Iso-pentane

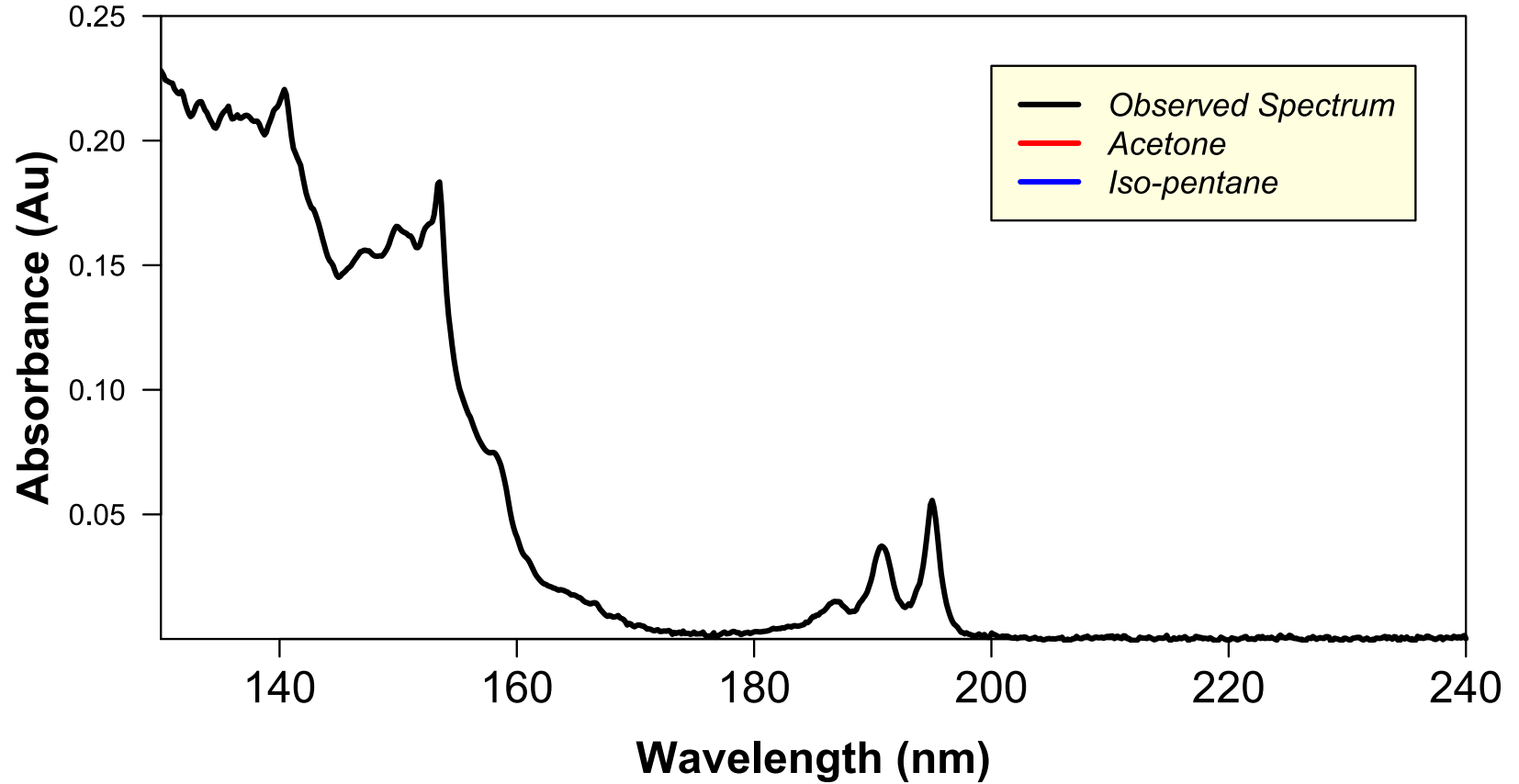

Coelution of Iso-pentane and Acetone

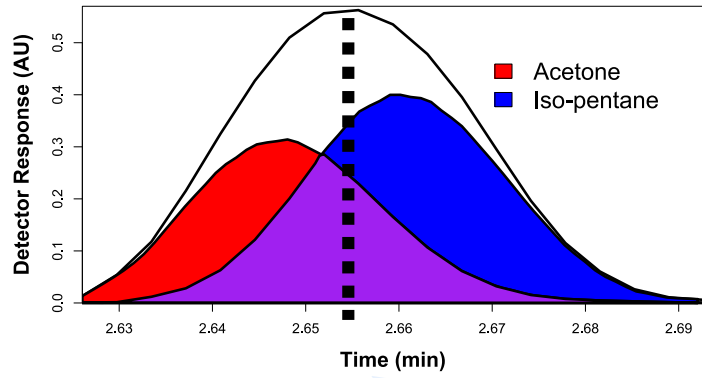

Spectral Absorbance of Acetone and Iso-pentane

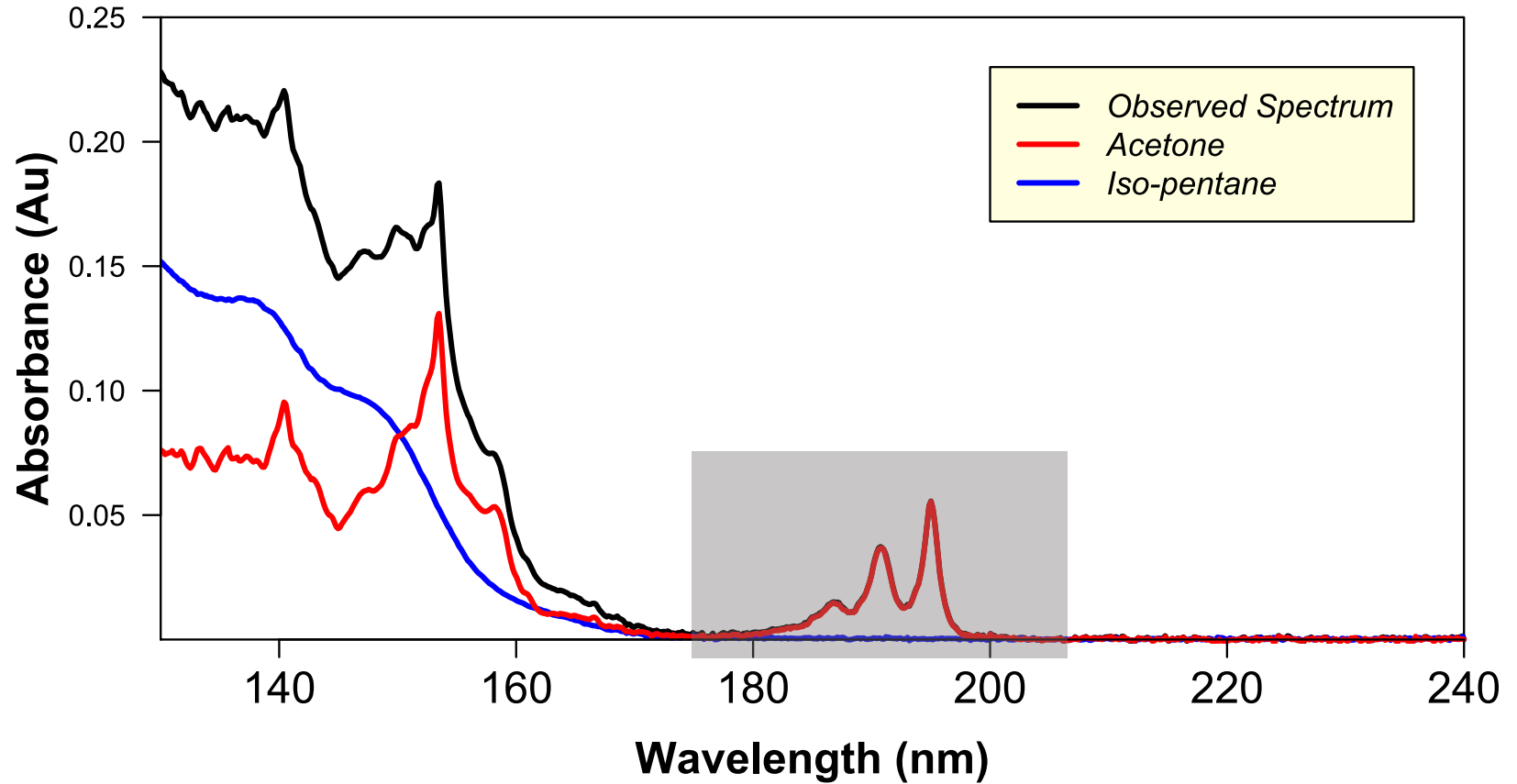



Designation: D8071 – 17

**Standard Test Method for
Determination of Hydrocarbon Group Types and Select
Hydrocarbon and Oxygenate Compounds in Automotive
Spark-Ignition Engine Fuel Using Gas Chromatography with
Vacuum Ultraviolet Absorption Spectroscopy Detection (GC-
VUV)¹**



✓ **Under 35 minute
run time**

✓ **Identification by
absorption and
retention time**

✓ **Spectral
deconvolution**

✓ **Easily add
compounds to
library**







Designation: D8071 – 17




Can we use ASTM D8071 conditions to analyze NTGAs?

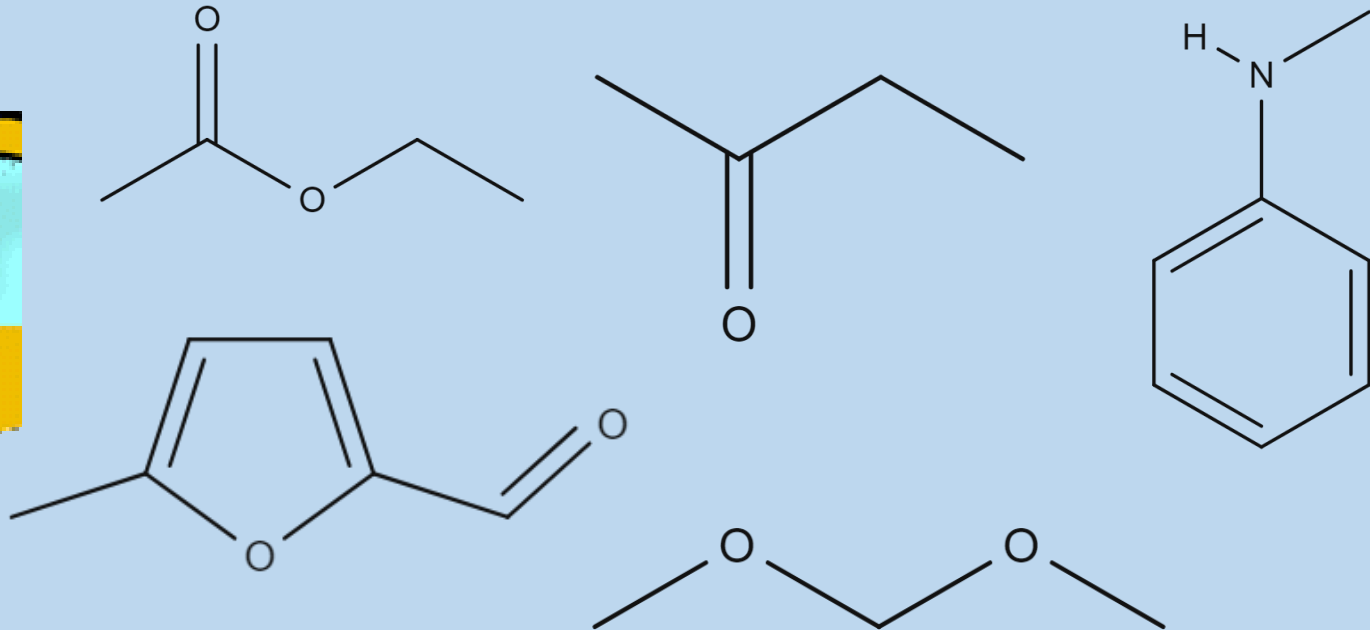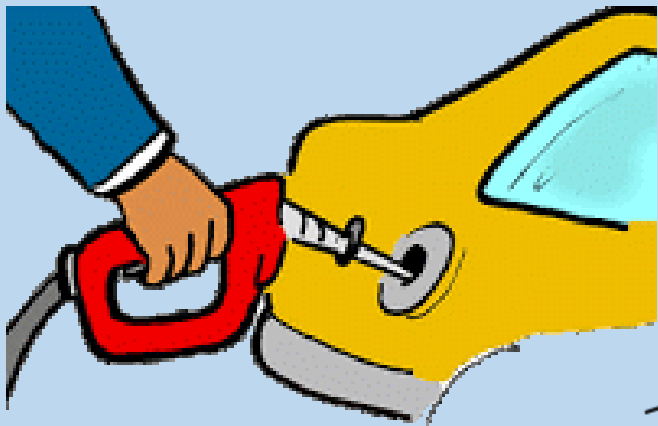

absorption and retention time

library



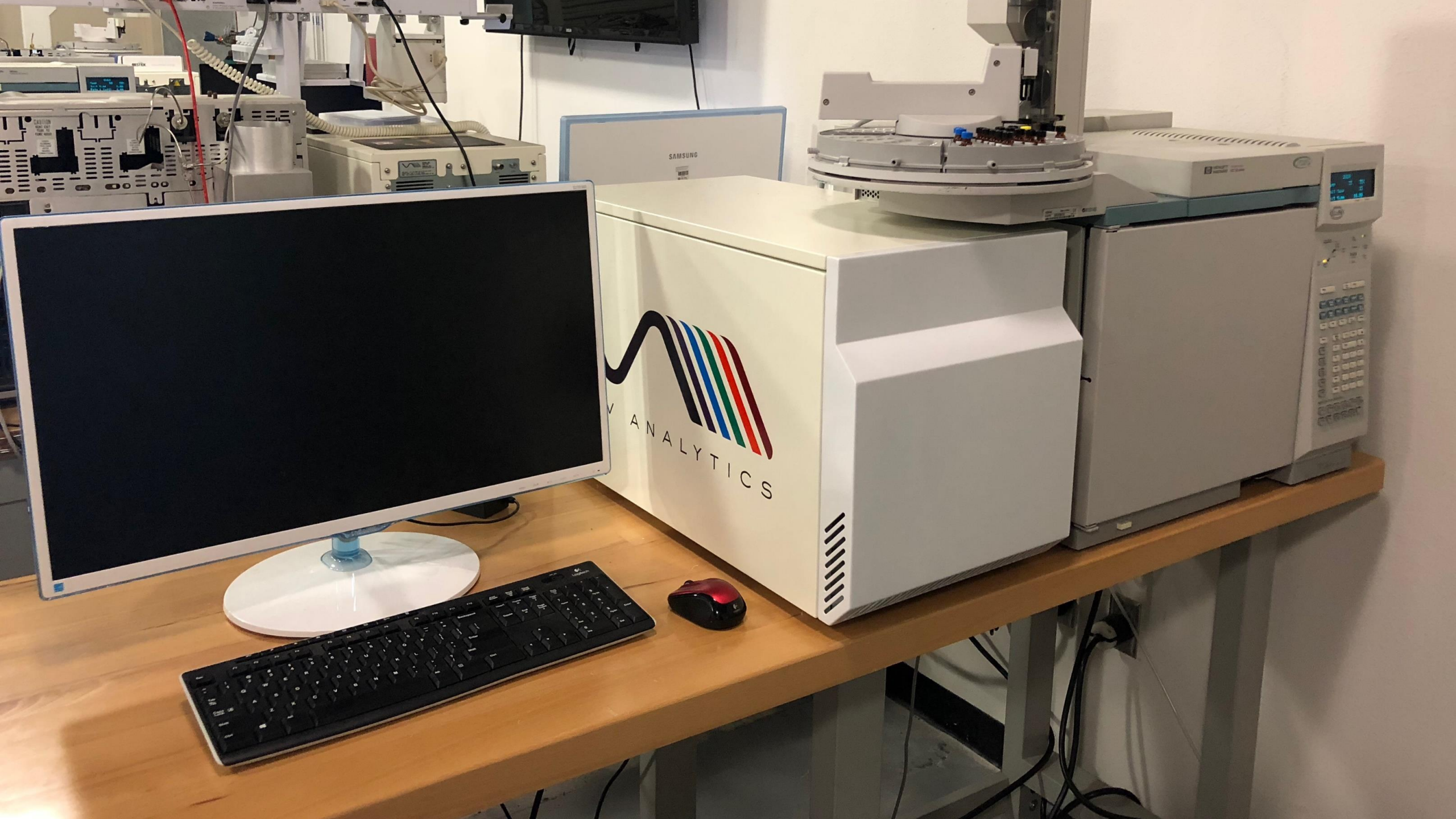

Experimental setup

1

2

3

Volumetric standards prepared in oxygenate-free gasoline

Experimental setup

1

Volumetric standards prepared in oxygenate-free gasoline

Acetone	Isobutyl acetate
Dimethoxymethane	3-Hexanone
Methyl acetate	Diethyl carbonate
2-Butanone	2,4-Dimethyl-3-pentanone
2-Methylfuran	Ethyl butanoate
Dimethyl carbonate	n-Butyl acetate
Ethyl acetate	Methyl pentanoate
2-Methyltetrahydrofuran	Ethyl pentanoate
2-Pentanone	Anisole
2-Pentanol	Isobutyl isobutyrate
2,5-Dimethylfuran	gamma-Valerolactone
4-Methyl-2-Pentanone	Aniline
2-Methyl-1-butanol	Methyl levulinate
sec-Butyl acetate	Ethyl levulinate
Cyclopentanone	N-Methylaniline

1%

3%

5%

10%

Volume %

Experimental setup

1

Volumetric standards prepared in oxygenate-free gasoline

Acetone	Isobutyl acetate
Dimethoxymethane	3-Hexanone
Methyl acetate	Diethyl carbonate
2-Butanone	2,4-Dimethyl-3-pentanone
2-Methylfuran	Ethyl butanoate
Dimethyl carbonate	n-Butyl acetate
Ethyl acetate	Methyl pentanoate
2-Methyltetrahydrofuran	Ethyl pentanoate
2-Pentanone	Anisole
2-Pentanol	Isobutyl isobutyrate
2,5-Dimethylfuran	gamma-Valerolactone
4-Methyl-2-Pentanone	Aniline
2-Methyl-1-butanol	Methyl levulinate
sec-Butyl acetate	Ethyl levulinate
Cyclopentanone	N-Methylaniline

Additional samples:

0.05%	0.1%	0.2%	0.5%	
1%	2%	5%	10%	20%

Experimental setup

1

2

3

**Analyzed with
ASTM D8071
using GC-VUV**

Experimental setup

ASTM D8071

Agilent 6890 GC

- Column: 30m x 0.25mm x 0.25 μ m Rxi-1ms
- GC Inlet: 250°C, split 300:1
- Constant Flow Mode: 1.0 mL/min He
- Injection volume: 1 μ L
- Oven: 35°C (10 min), 7°C/min to 200°C (0 min)

VUV Analytix VGA-100

- Makeup gas: 0.25 psi N₂
- Acquisition range: 125-240 nm
- Acquisition rate: 4.5 Hz
- Detector flow cell: 275°C
- Transfer line temperature: 275°C

Experimental setup

1



2



3

A = Absorption
 ϵ = Molar Absorptivity
b = Path Length
c = Concentration


Quantified

Beer-Lambert Law:

$$A = \epsilon bc$$

NTGA Compound	10%	5%	3%	1%	NTGA Compound	10%	5%	3%	1%
<i>Acetone</i>	10.9	4.96	3.00	0.956	<i>Isobutyl acetate</i>	10.5	5.40	3.20	1.04
<i>Dimethoxymethane</i>	10.4	5.12	3.13	1.08	<i>3-Hexanone</i>	10.0	5.25	3.25	1.09
<i>Methyl acetate</i>	10.4	5.20	3.14	1.06	<i>Diethyl carbonate</i>	10.5	5.36	3.18	0.954
<i>2-Butanone</i>	10.4	5.12	3.09	1.11	<i>2,4-Dimethyl-3-pentanone</i>	10.1	5.15	3.18	1.08
<i>2-Methylfuran</i>	10.2	5.24	3.17	1.05	<i>Ethyl butanoate</i>	10.0	5.22	3.19	1.06
<i>Dimethyl carbonate</i>	10.5	5.21	3.09	1.05	<i>n-Butyl acetate</i>	10.2	5.27	3.16	1.06
<i>Ethyl acetate</i>	10.3	5.65	3.66	1.25	<i>Methyl pentanoate</i>	10.1	5.17	3.13	1.03
<i>2-Methyltetrahydrofuran</i>	10.2	5.13	3.13	1.04	<i>Ethyl pentanoate</i>	10.3	5.37	3.31	1.18
<i>2-Pentanone</i>	10.6	5.32	3.18	1.01	<i>Anisole</i>	10.1	5.16	3.03	1.04
<i>2-Pentanol</i>	10.2	5.19	3.32	1.15	<i>Isobutyl isobutyrate</i>	10.0	5.12	3.13	1.04
<i>2,5-Dimethylfuran</i>	10.2	5.18	3.18	1.03	<i>gamma-Valerolactone</i>	6.71	5.02	3.08	1.00
<i>4-Methyl-2-Pentanone</i>	10.3	5.30	3.19	1.06	<i>Aniline</i>	10.3	5.23	3.19	1.07
<i>2-Methyl-1-butanol</i>	10.3	5.38	3.21	1.08	<i>Methyl levulinate</i>	10.2	5.13	3.19	1.10
<i>sec-Butyl acetate</i>	10.5	5.43	3.25	1.08	<i>Ethyl levulinate</i>	10.2	5.20	3.16	1.06
<i>Cyclopentanone</i>	10.1	5.18	3.14	1.00	<i>N-Methylaniline</i>	10.3	5.32	3.28	1.08

NTGA Compound	10%	5%	3%	1%	NTGA Compound	10%	5%	3%	1%
Acetone	10.9	4.96	3.00	0.956	Isobutyl acetate	10.5	5.40	3.20	1.04
Dimethoxymethane	10.13	5.13	3.13	1.08	3-Hexanone	10.0	5.25	3.25	1.09
Methyl acetate	10.6	5.06	3.06	0.96	Diethyl carbonate	10.5	5.36	3.18	0.954
2-Butanol	10.1	5.1	3.1	1.0	2,4-Dimethyl-3-pentanone	10.1	5.15	3.18	1.08
2-Methyl-2-butanol	10.1	5.1	3.1	1.0	Ethyl butanoate	10.0	5.22	3.19	1.06
Dimethyl ether	10.1	5.1	3.1	1.0	n-Butyl acetate	10.2	5.27	3.16	1.06
Ethyl acetate	10.1	5.1	3.1	1.0	Methyl pentanoate	10.1	5.17	3.13	1.03
2-Methyltetrahydrofuran	10.1	5.1	3.1	1.0	Ethyl pentanoate	10.3	5.37	3.31	1.18
2-Pentanone	10.1	5.1	3.1	1.0	Anisole	10.1	5.16	3.03	1.04
2-Pentanol	10.2	5.19	3.32	1.15	Isobutyl isobutyrate	10.0	5.12	3.13	1.04
2,5-Dimethylfuran	10.2	5.18	3.18	1.03	gamma-Valerolactone	6.71	5.02	3.08	1.00
4-Methyl-2-Pentanone	10.3	5.30	3.19	1.06	Aniline	10.3	5.23	3.1	1.07
2-Methyl-1-butanol	10.3	5.38	3.21	1.08	Methyl levulinate	10.2	5.13	3.19	1.10
sec-Butyl acetate	10.5	5.43	3.25	1.08	Ethyl levulinate	10.2	5.20	3.16	1.06
Cyclopentanone	10.1	5.18	3.14	1.00	N-Methylaniline	10.3	5.32	3.28	1.08

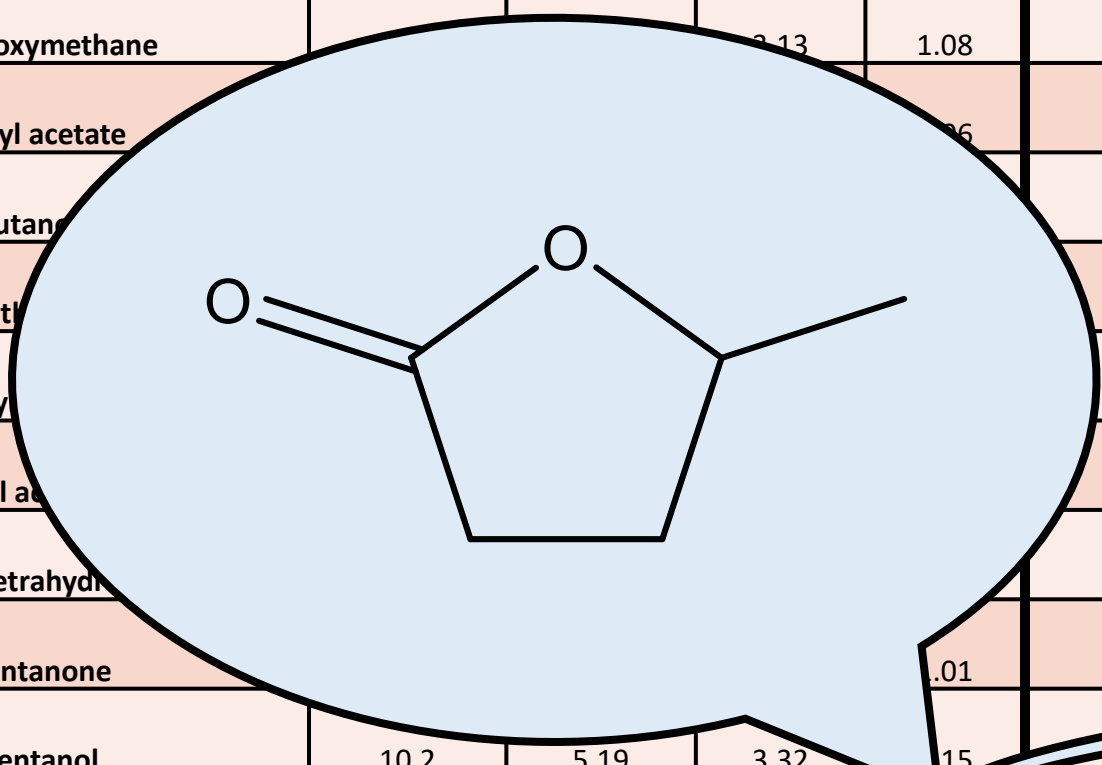


gamma-Valerolactone 6.71 5.02 3.08 1.00


Oxygenate-Free Gasoline


NTGA Compound	10%	5%	3%	1%
gamma-Valerolactone	<u>6.71</u>	5.02	3.08	1.00

gamma-Valerolactone

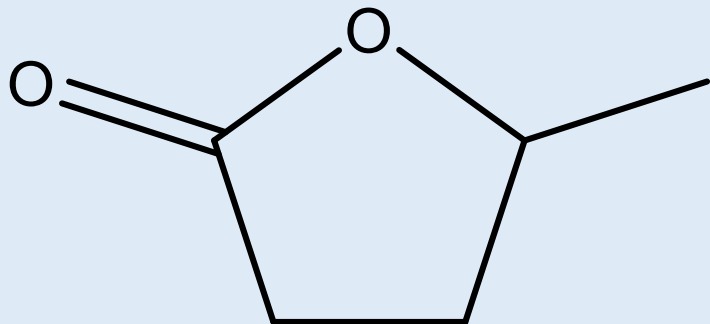

Insoluble

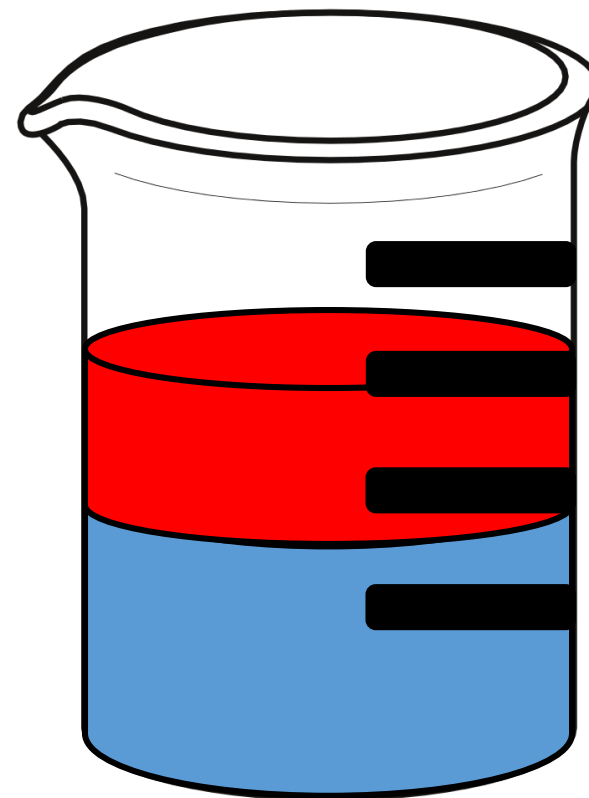

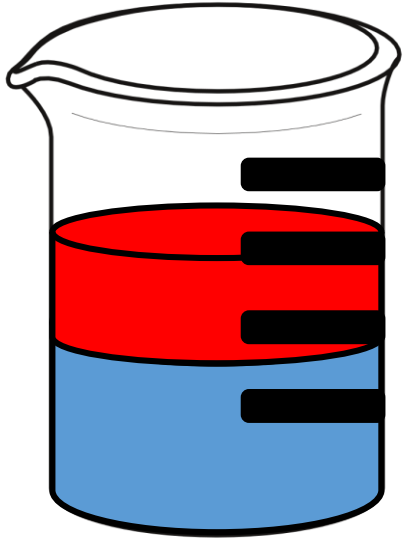


(No ethanol)
Insoluble



(10% ethanol)
Soluble


Oxygenate-Free Gasoline				
NTGA Compound	10%	5%	3%	1%
gamma-Valerolactone	<u>6.71</u>	5.02	3.08	1.00

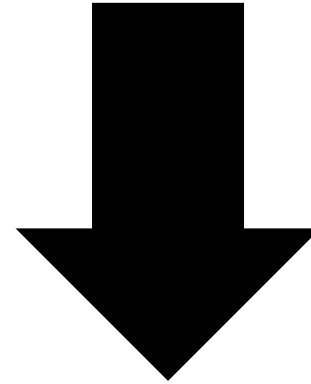


Murphy Express Gasoline				
NTGA Compound	10%	5%	3%	1%
gamma-Valerolactone	<u>10.3</u>	5.51	3.26	1.05

NTGA Compound	20%	10%	5%	2%	1%	0.5%	0.2%	0.1%	0.05%
<i>Dimethoxymethane</i>	20.9	9.85	4.90	1.83	1.00	0.49	0.22	0.09	0.06
<i>N-Methylaniline</i>	22.8	11.2	5.62	2.20	1.11	0.54	0.22	0.07	0.03
<i>2,5-Dimethylfuran</i>	20.7	9.96	4.88	1.83	1.01	0.47	0.16	0.06	0.02


Dimethoxymethane

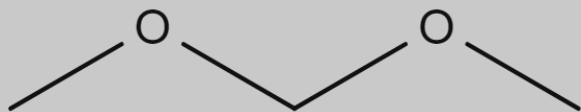


N-Methylaniline

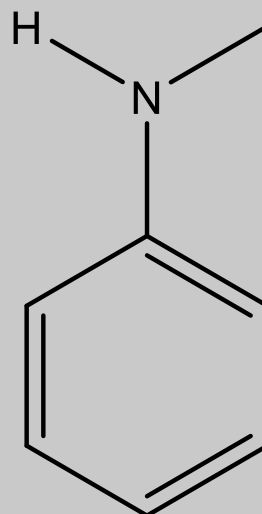


2,5-Dimethylfuran

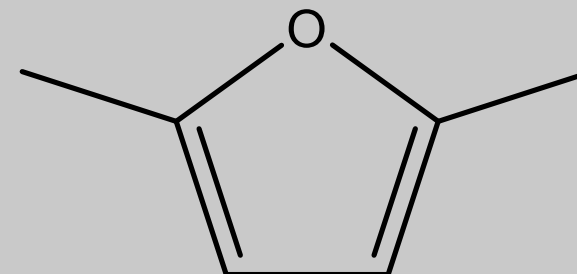


Dimethoxymethane in Gasoline



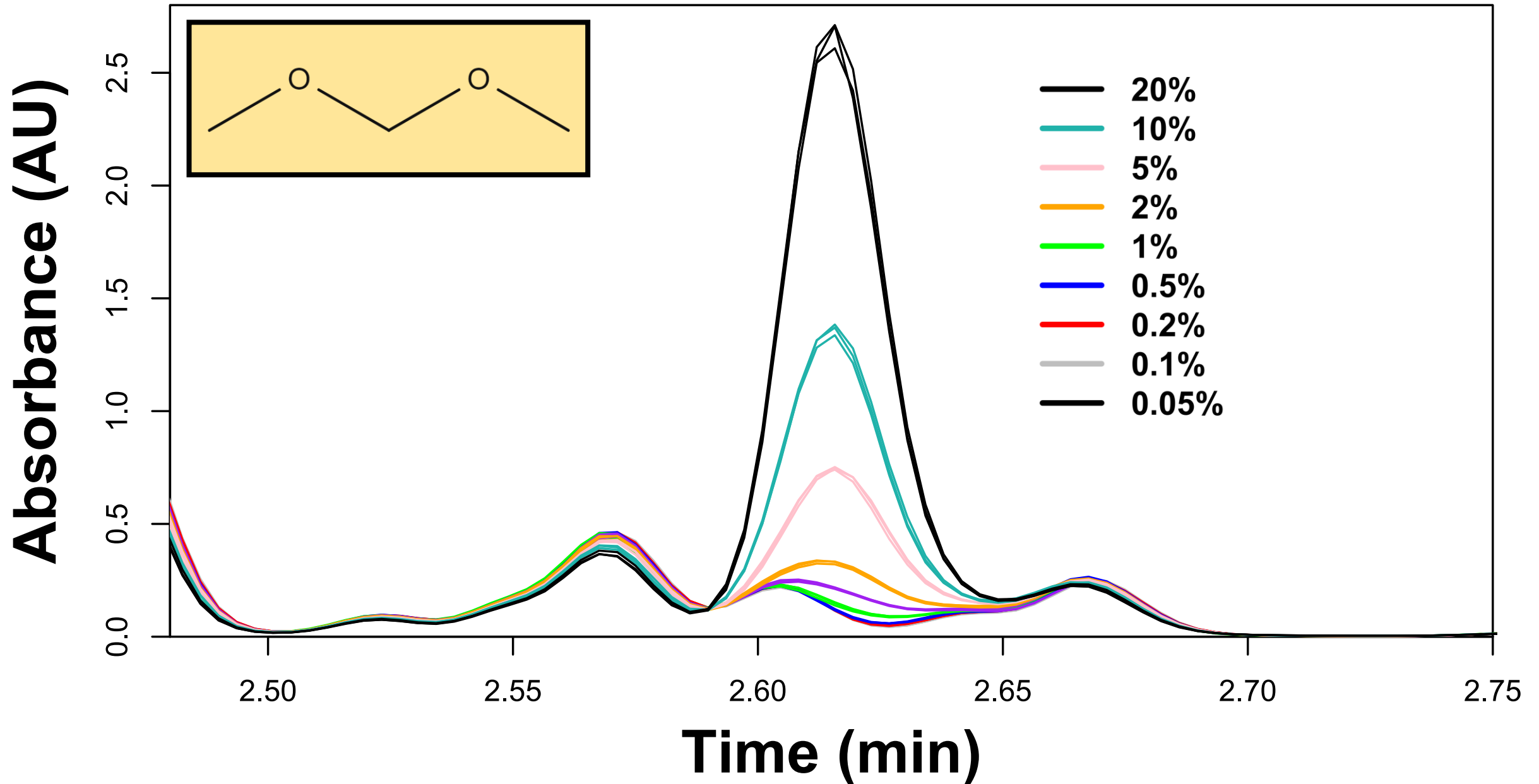





Dimethoxymethane Response vs. Concentration %

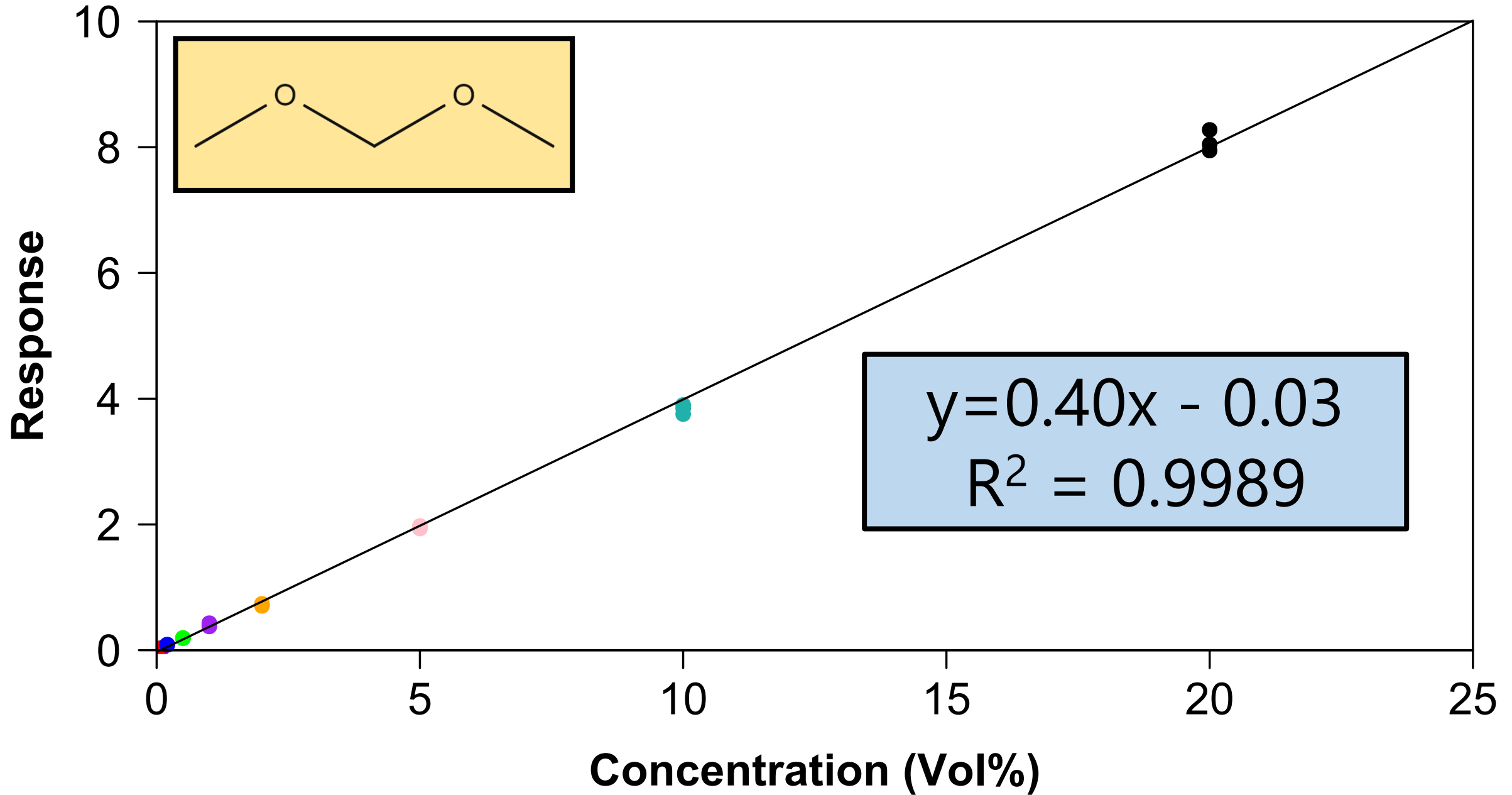


10% Ethyl Acetate in Gasoline

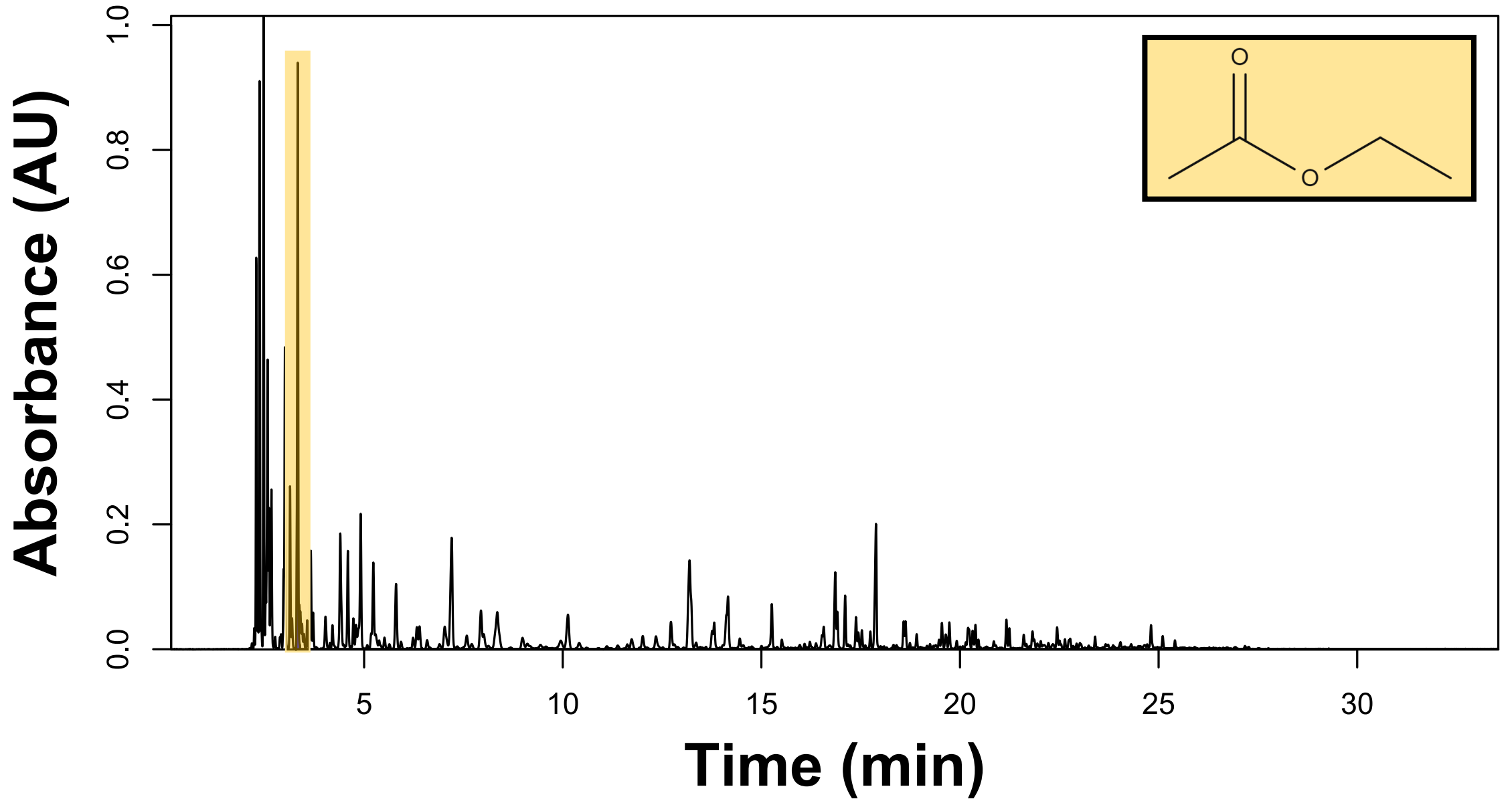

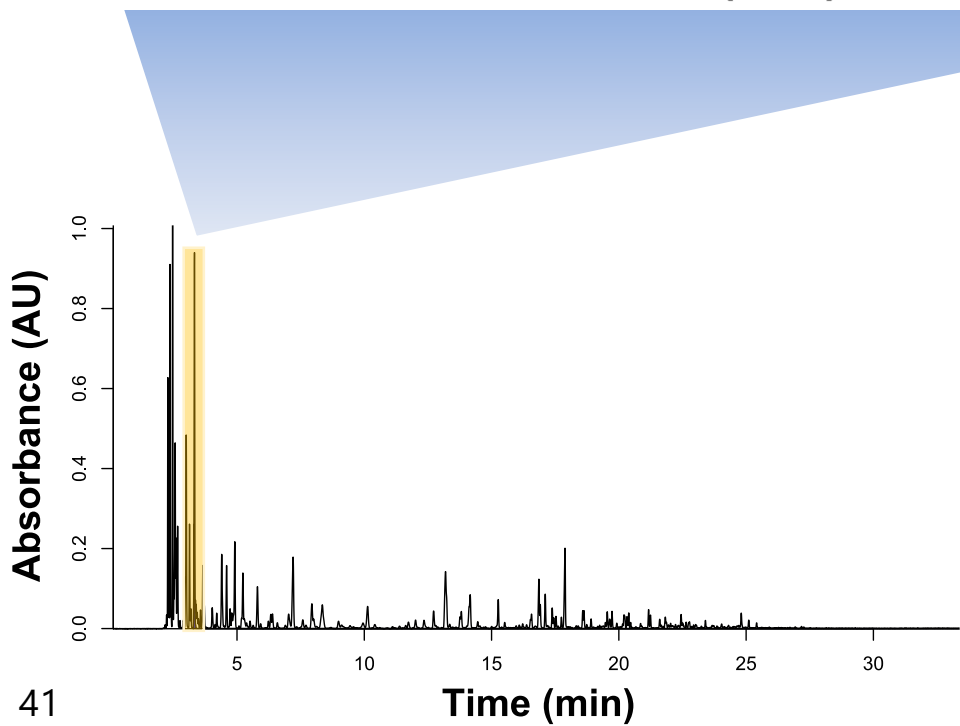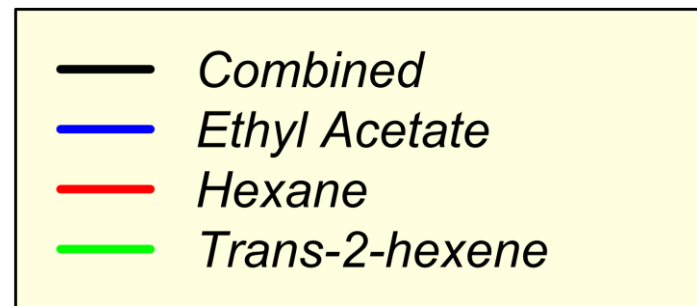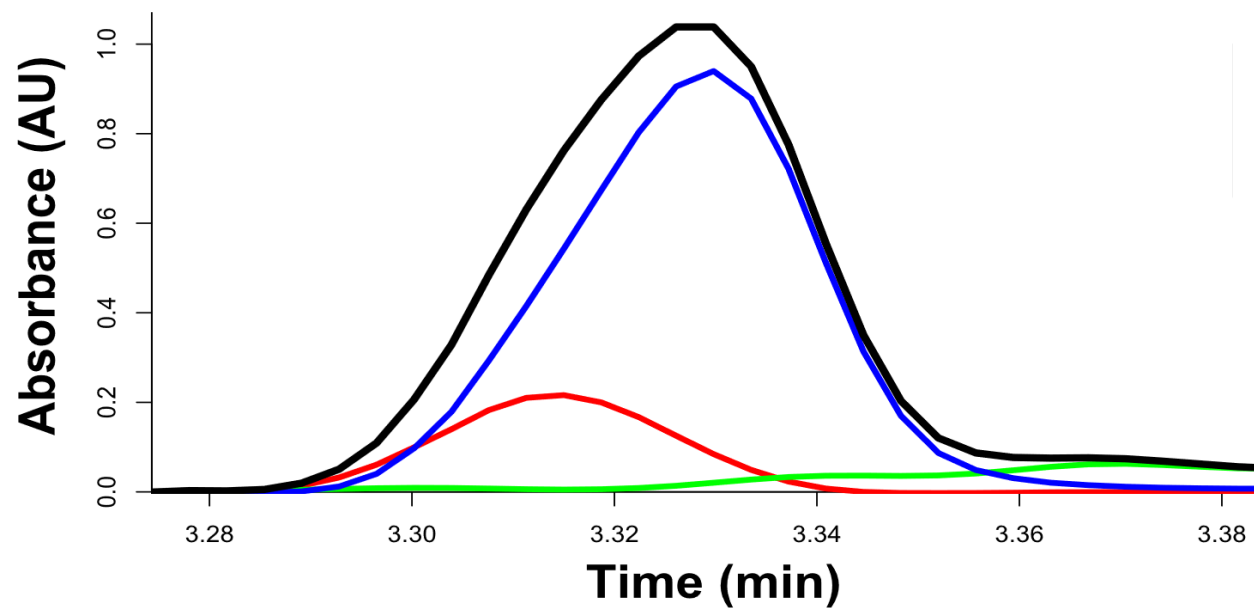

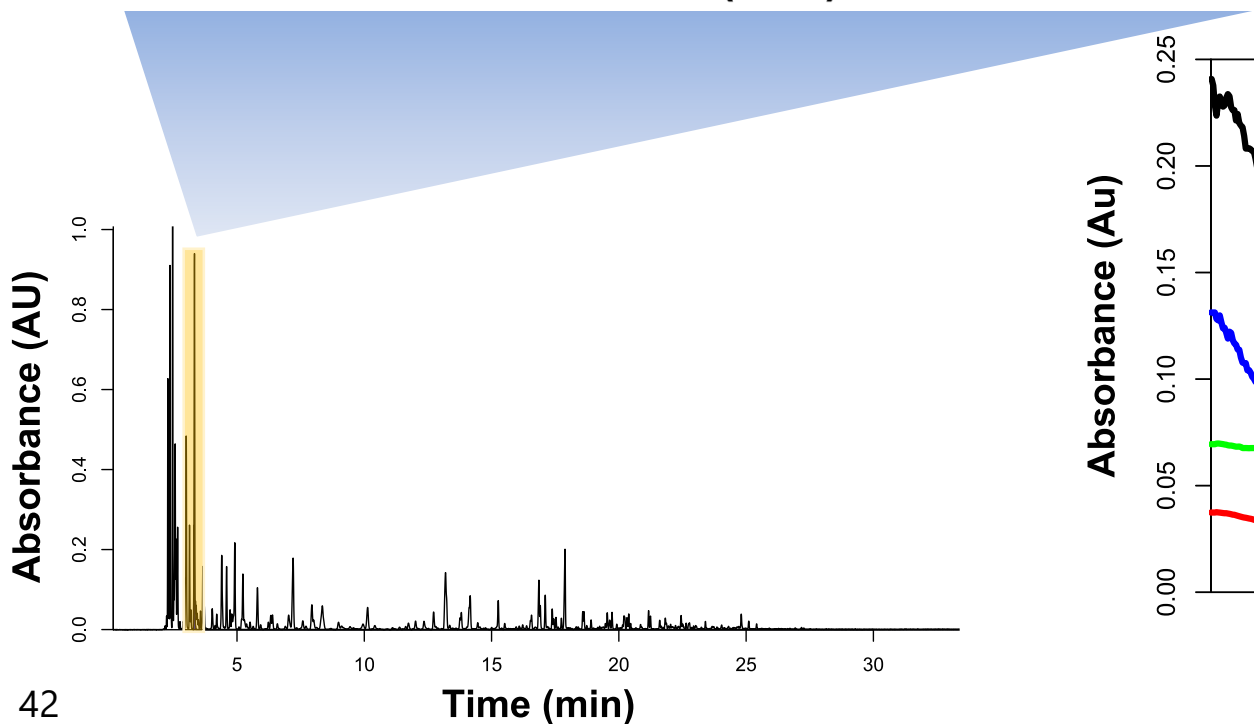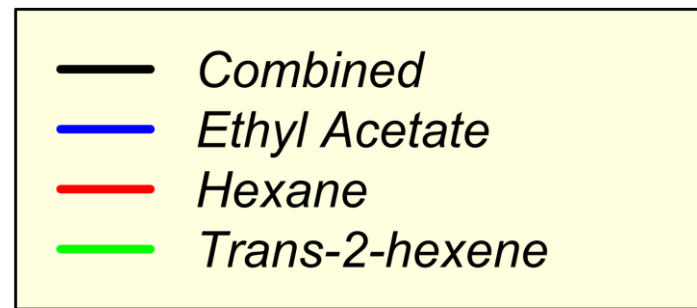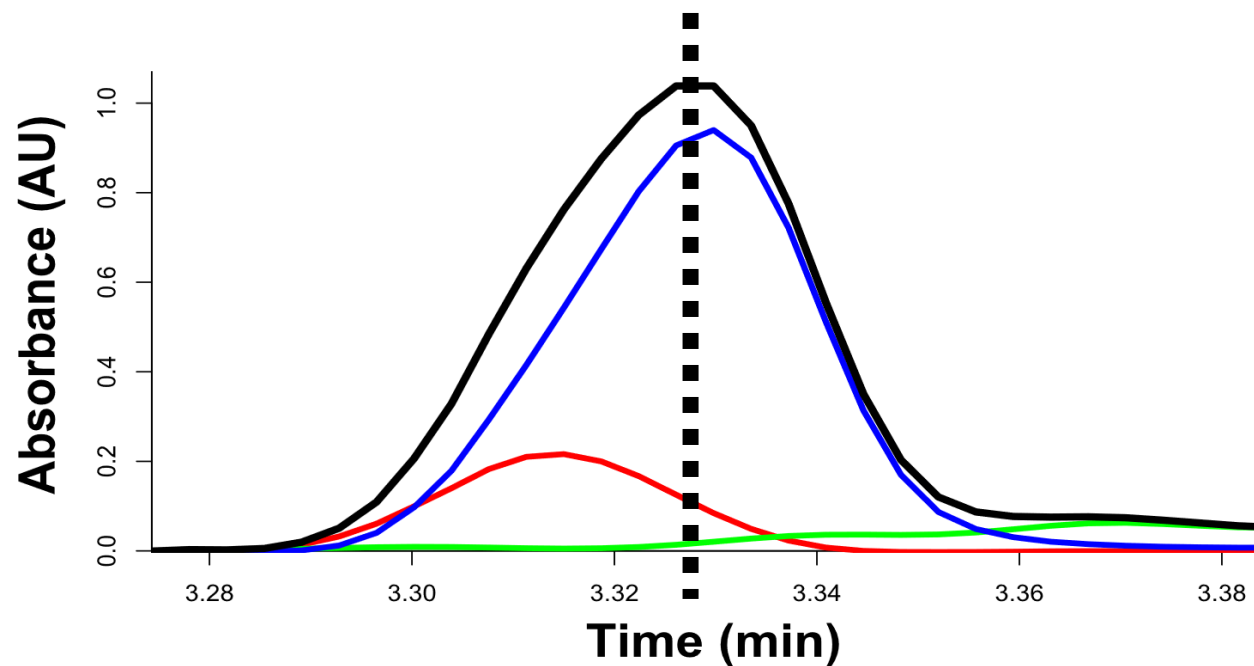


Spectral Deconvolution of Ethyl Acetate

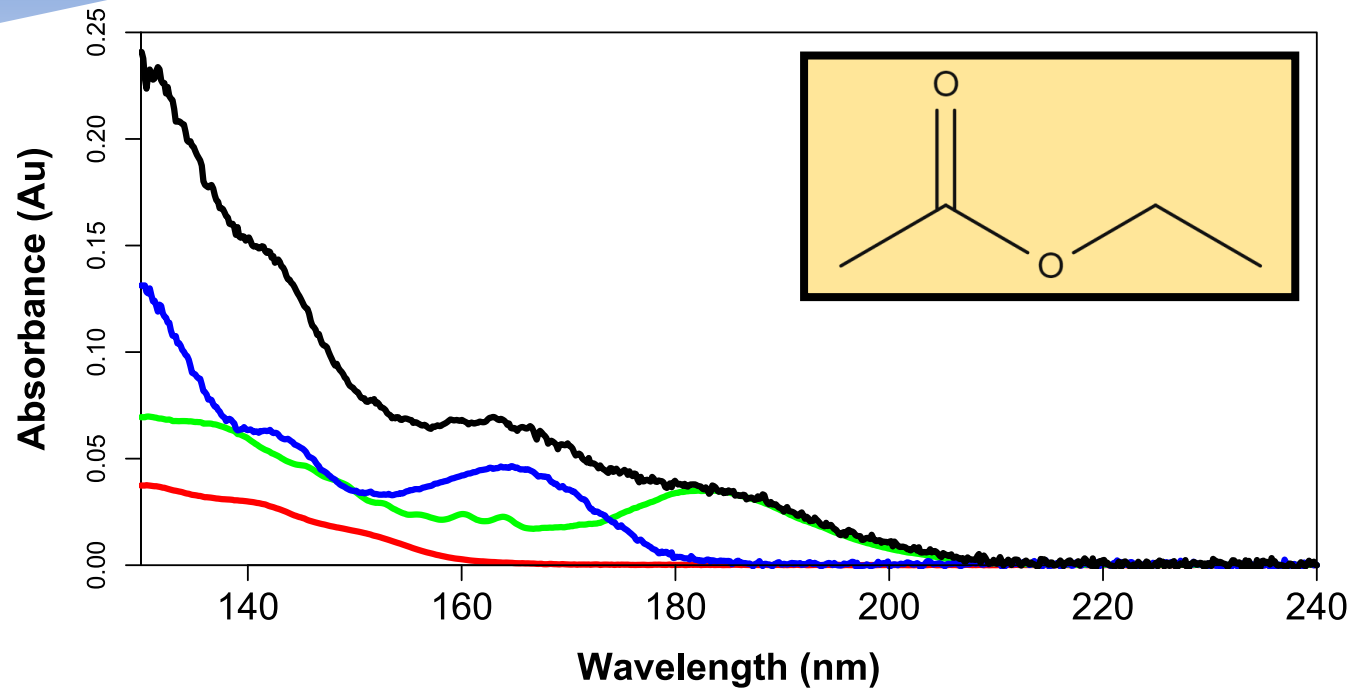

Conclusions

- NTGAs can be analyzed using GC-VUV and ASTM D8071 conditions
 - Compounds easily added to library
 - Deconvolve coeluting compounds
 - Accurately quantified
- ASTM D8071 conditions are favorable over ASTM D6730
 - Shorter run time
 - Confidence in identification
 - No method changes necessary

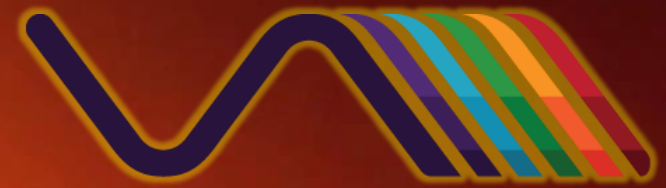

VUV ANALYTICS



Image:
<https://www.videoblocks.com/video/oil-platform-at-sea-at-sunset-lkqbl2d>

Questions?

Ryan.Schonert@VUVanalytics.com